

Application Data Sheet

No.101

GC-MS Gas Chromatograph Mass Spectromete

Analysis of Toxicological Substances in Whole Blood Using the Smart Forensic Database (2)

Smart Forensic Database is an MRM database containing retention indices, MRM transitions, collision energies, and quantitation/confirmation ion ratios for 201 toxicological substances often involved in poisonings. Smart MRM, which is provided with the GCMS-TQ8040, can create analysis methods for Scan/MRM and MRM automatically from the database.

This article introduces an example of using a simultaneous Scan/MRM measurement method created from the Smart Forensic Database to measure a whole-blood sample, and then applying the GC/MS Forensic Toxicology Database to the scan data obtained.

GC/MS Forensic Toxicological Database

The GC/MS Forensic Toxicological Database is intended for Scan analysis, and is registered with 726 toxicological compounds (1476 types). The database contains toxicological substance names, quantitation and confirmation ions, retention indices, and standard mass spectra. The retention times for the registered toxicological substances are accurately estimated simultaneously from low-boiling point components to high-boiling point components, using measurement data from a standard n-alkane mixture via the GCMSsolution AART function. It is easy to determine whether samples contain toxicological substances by using the mass chromatograms for the quantitation and confirmation ions contained in the database.

Experimental

Liquid-liquid extraction via EXtrelut[®] NT3 was used to pretreat the whole-blood sample. The collected wholeblood sample was measured into 1 mL portions for acidic fractionation and basic fractionation, and each portion was diluted with 1 mL of Milli-Q water. The acidic fraction was adjusted to a pH 5 using 10 % hydrochloric acid, and the basic fraction was adjusted to a pH 9 using 10 % ammonia water. The respective solutions were added to the EXtrelut[®] NT3 columns and left to stand for 30 minutes, after which each was eluted with a 10 mL chloroform:isopropanol (3:1) mixture. The extracted solutions of acidic fraction and basic fraction was re-dissolved in a 200 μ L chloroform:isopropanol (3:1) mixture.

Analytical Conditions

The conditions registered in Smart Forensic Database were used as the GC-MS/MS analysis conditions. For the compounds subject to MRM measurement, a simultaneous Scan/MRM analysis method was created, in which the 201 components registered in the database were set.

Table 1: Analytical Conditions

GC-MS: GCMS-TQ8040 Column: Rxi®-5SilMS (Length 30 m, 0.25 mm l.D., df=0.25 μm) Glass liner: Splitless insert with wool (PN: 221-48876-03)					
[GC] Injection temp.: Column oven temp.: Injection mode: Flow control mode: Injection volume:	260 °C 60 °C (2 min) \rightarrow (10 °C /min) \rightarrow 320 °C (15 min) Splitless Linear velocity (45.6 cm/sec) 1 µL	[MS] Interface temp.: Ion source temp.: Acquisition mode: Scan event time: Scan mass range: Scan speed: MRM event time: Total loop time:	280 °C 200 °C Scan/MRM 0.1 sec m/z 43 – 600 10,000 u/sec 0.3 sec 0.4 sec		

Results

The extracted whole-blood samples were measured in the simultaneous Scan/MRM analysis mode. By analyzing the scan data using the GC/MS Forensic Toxicological Database, pentedrone, a type of cathinone, was identified (Fig. 1). If a simultaneous Scan/MRM analysis is done, the obtained MRM data can be analyzed via Smart Forensic Database to identify toxicological substances often involved in poisonings, while the scan data can be analyzed to confirm identifications using the mass spectra, or to screen for approximately 1400 toxicological substances using the GC/MS Forensic Toxicological Database.

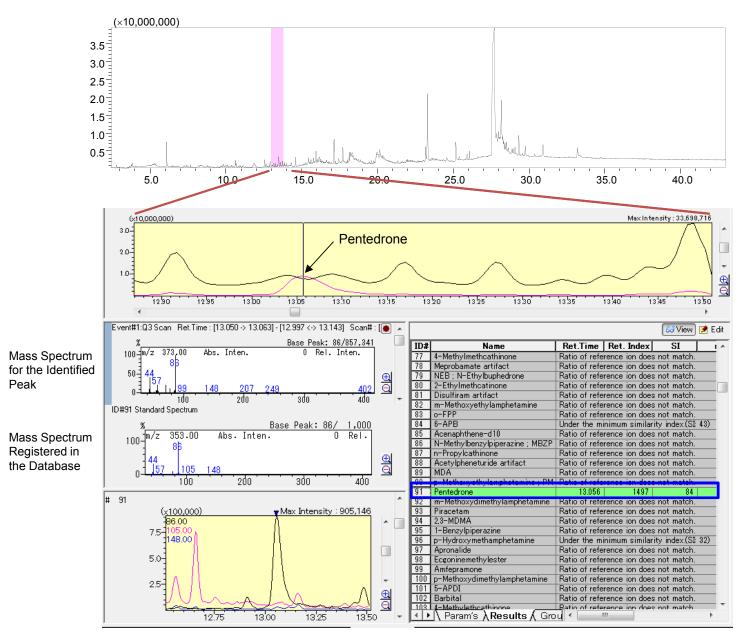


Fig. 1: Results from an Analysis of the Scan Data from a Whole-Blood Sample Using the GC/MS Forensic Toxicological Database

This data was provided by Associate Professor Kei Zaitsu in the Department of Legal Medicine & Bioethics, Nagoya University Graduate School of Medicine.

First Edition: September, 2014



Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.