

Application
Data Sheet

No. 63

System Gas Chromatograph

**Trace H₂ Analysis System
Nexis GC-2030TH2
GC-2014TH2**

The system enables a quantitative and qualitative analysis of H₂ in municipal gas. A total of 1 valve and 2 columns are used in this GC system. N₂ is used as the carrier gas. Sample is introduced into one sample loop for determination. Using a pre-column, C₂-C₃ components are back-flushed. The valve timing then allows O₂ and N₂ to directed to molecular sieve column for separation and TCD for detection. The analysis time is approximately 4 minutes. LabSolutions chromatography workstation system handles all aspects of GC control, automation, and data handling.

Analyzer Information

System Configuration:

One valve / Two packed columns with one TCD detector

Sample Information:

H₂

Methods met:

ASTM-D2504

Concentration Range:

No.	Name of Compound	Concentration Range		Detector
		Low Conc.	High Conc.	
1	H ₂	5ppm	500ppm	TCD-1

Detection limits may vary depending on the sample. Please contact us for more consultation.

System Features

- 5 minutes analysis for H₂ analysis can be carried out
- Single TCD channel
- Good repeatability

Typical Chromatograms

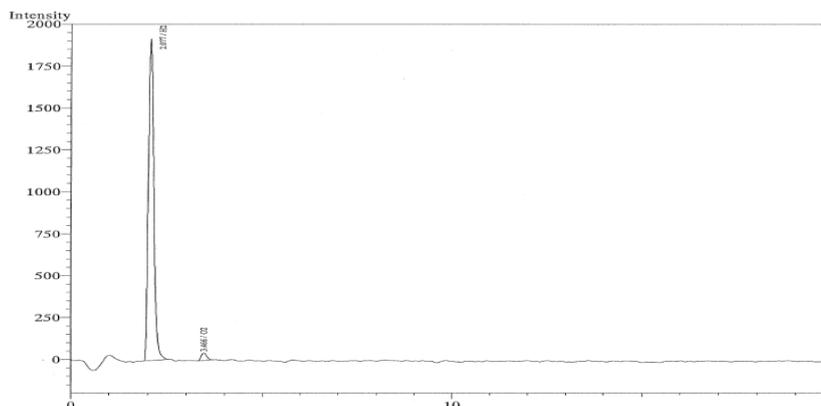


Fig. Chromatogram of TCD

First Edition: November, 2017