

Application News

Liquid Chromatograph Mass Spectrometry

No. **C166**

Analysis of Residual Pesticides in Vegetable Extracts Using LC/MS/MS <LCMS-8050>

In order to prevent residual pesticides and chemicals in food items from adversely affecting human health, the maximum residue limit of all pesticides, feed additives, and veterinary pharmaceuticals is set for each food item within the range of amounts that were determined to be safe for people to consume (the Positive List System). The Food Sanitation Act prohibits selling and importing food items with amounts of pesticides or other chemicals that exceed the maximum residue limit. Furthermore, maximum usage limits of pesticides and other chemicals are prescribed in the Agricultural Chemicals Regulation Law to prevent them from remaining in food items at amounts that exceed the limits.

This article introduces an analysis of pesticides contained in green onion extracts using liquid chromatography tandem mass spectrometry (LC/MS/MS). The analyzed pesticides include endosulfan (benzoepin) which is a substance that was added on April 1, 2012 to the Ministerial Ordinance on Prohibition of Agricultural Chemical Sales.

While endosulfan has conventionally been analyzed by gas chromatography mass spectrometry (GC/MS), detection by liquid chromatography mass spectrometry (LC/MS) with high sensitivity has become possible by optimizing the temperature and other conditions of the electrospray ionization interface for LC/MS. Good recovery factor resulted even for the green onion extracts obtained by the QuEChERS method which is known to have a relatively strong matrix influence. The example analyses introduced here cover endosulfan, thiamethoxam, dinotefuran, and an iprodione metabolite in green onion extract as well as imazalil, fludioxonil, pyrimethanil, and azoxystrobin in orange extract. M. Tanaka

Analytical Conditions

The analytical conditions of high performance liquid chromatography (HPLC) and Mass Spectrometry (MS) are shown in Tables 1 and 2.

Table 1 Analytical Conditions for HPLC

| Column | : Shim-pack [™] XR-ODS II |
|------------------|---|
| | (75 mm L.×2.1 mm l.D., 2.2 μm) |
| Mobile phase | : A) 0.5 mmol/L ammonium acetate-water |
| | B) 0.5 mmol/L ammonium acetate-methanol |
| Time programs | : B conc. 10 % (0 min) $ ightarrow$ 95 % (4.50-8.00 min) $ ightarrow$ |
| | 10 % (8.51-12.00 min) |
| Column temp. | : 40 °C |
| Flow rate | : 0.2 mL/min |
| Injection Volume | : Green Onion 1 μL, Orange 2 μL |
| | |

Table 2 Analytical Conditions for MS

| | Green Onion | Orange | |
|---------------------|-------------------------|-------------------------|--|
| Probe Voltage | 0.5 kV (ESI-Positive) / | 0.5 kV (ESI-Positive) / | |
| | –0.5 kV (ESI-Negative) | –0.5 kV (ESI-Negative) | |
| DL Temp. | 150 °C | 150 °C | |
| Block Heater Temp. | 300 °C | 500 °C | |
| Interface Temp. | 100 °C | 400 °C | |
| Nebulizing gas flow | 3.0 L/min | 3.0 L/min | |
| Drying gas flow | 13.0 L/min | 10.0 L/min | |
| Heating gas flow | 7.0 L/min | 10.0 L/min | |

MRM Transition Conditions

Data of the multiple reaction monitoring (MRM) transition and the calibration curve of each pesticide are shown in Table 3. An acceptable coefficient of determination was obtained for each pesticide in the calibration curve concentration range with each value being higher than 0.99.

| Pesticide | Polarity (+/–) | Precursor (<i>m/z</i>) | Product (<i>m/z</i>) | Calibration Curve Conc. Range (ng/mL) | Coefficient of Determination (r ²) |
|-----------------------------------|-------------------|-----------------------------|---------------------------|---|--|
| Dinotefuran | + | 203.10 | 129.2 | 0.05-100 | 0.9981 |
| Thiamethoxam | + | 292.10 | 211.0 | 0.05-100 | 0.9978 |
| Dinotefuran | - | 201.00 | 61.0 | 0.5-100 | 0.9979 |
| Iprodione | - | 328.05 | 141.0 | 0.02-100 | 0.9992 |
| metabolite | | | | | |
| Endosulfan (α + β) | - | 404.70 | 268.9 | 0.2-100 | 0.9996 |
| Azoxystrobin | + | 404.10 | 372.1 | 0.01-50 | 0.9999 |
| Pyrimethanil | + | 200.20 | 107.1 | 0.02-50 | 0.9981 |
| Imazalil | + | 296.90 | 159.0 | 0.05-50 | 0.9949 |
| Fludioxonil | - | 246.90 | 180.1 | 0.05-50 | 0.9995 |

Linearity of Calibration Curves

The calibration curves of common pesticides, an iprodione metabolite and endosulfan (α + β), are shown in Figs. 1 and 2.



Fig. 1 Calibration Curve (Iprodione Metabolite)



Fig. 2 Calibration Curve (Endosulfan $(\alpha+\beta)$)

Chromatograms of Samples

The chromatograms of samples are shown in Figs. 3 and 4. Through selective analysis in which impurity peaks are not

detected, each pesticide was well separated and chromatograms with good peak shapes were obtained.









Recovery Factor

Recovery factor of all pesticides in the green onion and orange extracts are shown in Table 4.

The recovery factor was obtained using the following formula:

Recovery factor (%) = ((Measured concentration – Concentration of blank sample) / Actual spike concentration) × 100 Good recovery factors were obtained for the green onion and orange extracts in the range of 88 to 104 % and 74 to 98 %, respectively.

The peak area repeatabilities (%RSD, n = 3) of the samples are shown in Table 4. The %RSD for each pesticide is favorable with all values being lower than 8 %.

| Pesticide | Spike Concentration (ng/mL) | Recovery Factor (%) | %RSD | Pesticide | Spike Concentration (ng/mL) | Recovery Factor (%) | %RSD |
|---------------------------------|--------------------------------|------------------------|------|--------------|--------------------------------|------------------------|------|
| Dinotefuran | 100 | 101 | 0.70 | Azoxystrobin | 2 | 93 | 2.66 |
| Thiamethoxam | 10 | 104 | 5.50 | Pyrimethanil | 40 | 74 | 5.08 |
| Iprodione metabolite | 25 | 88 | 2.80 | Imazalil | 10 | 98 | 0.66 |
| Endosulfan ($\alpha + \beta$) | 50 | 104 | 5.36 | Fludioxonil | 20 | 83 | 7.78 |

Table 4 Recovery Factor and Peak Area Repeatability

First Edition: May 2018



Shimadzu Corporation www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedure.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. Shimadzu disclaims any proprietary interest in trademarks and trade names used in this publication other than its own. See http://www.shimadzu.com/about/trademarks/index.html for details.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.