

Liquid Chromatography Mass Spectrometry

## Application News

## No.**C107**

Vitamin D is a fat-soluble vitamin that functions as a hormone in the body to regulate calcium metabolism. In addition, there is a report that vitamin D has a number of other physiological roles, including maintaining muscle strength, modulating immune function, and regulating blood pressure. Moreover, vitamin D functions to help regulate cellular differentiation, the biological process by which cells become specialized for a specific function.



Fig. 1 Structure of Vitamin D Metabolites

Quantitative Analysis of Vitamin D Metabolite Using Triple Quadrupole LC/MS/MS

> Due to these physiological functions, vitamin D has been implicated in protection against muscle pain and weakness, certain autoimmune diseases, hypertension, and even some forms of cancer.

> This sheet describes analysis of 25 hydroxy vitamin  $D_2$  and  $D_3$  in human serum using LC/MS/MS with ISOLUTE® SLE+.

[Sample Preparation Procedure]

1. Sample Pre-treatment:

To a mixture of 150  $\mu L$  water and 150  $\mu L$  isopropanol with ISTD add 100  $\mu L$  of sample and mix for 10 seconds.

2. Sample Load:

Load pre-treated sample (400  $\mu$ L) to plate followed by a pulse of positive pressure (2-3 seconds) to initiate flow and leave to flow under gravity for five minutes.

3. Analyte Elution:

Elute with heptane (700  $\mu$ L × 2) directly in to a deep well collection plate. Leave each aliquot to flow under gravity for 5 minutes then apply positive pressure to completely remove the final volume.

4. Post extraction:

Transfer extraction solvent to glass vials and evaporate to dryness at 40 °C.

Reconstitute in 100  $\mu$ L of solution containing: 33 % of mobile phase A and 67 % of mobile phase B v/v. Vortex the aliquot, centrifuge and transfer supernatant to new deep well plate.

[LC] NexeraX2 Syste Column Column Temp.	em :YMC Triart C18 (50 mm L. × 2 mm I.D., 1.9 μm) :40 °C
Mobile Phase A	Water : Methanol : Formic Acid = 50 : 50 : 0.025
Mobile Phase B	: Methanol
Time Program	:67 %B (0 - 1 min) - 67 %B - 95 %B (1 - 3 min) - 95 %B (3 - 4 min) - 67 %B (4 - 6 min)
Flowrate	: 0.5 mL/min
Injection Volume	: 10 µL
[MS] LCMS-8050	
Ionization	: ESI Positive
DL Temp.	: 150 ℃
Block Heater Temp.	: 400 °C
Interface Temp.	: 180 °C
Nebulizing Gas Flow	v : 3 L/min
Drying Gas Flow	: 10 L/min
Heating Gas Flow	:10 L/min

A seven point calibration curve was generated. The calibrant levels for the 25-hydroxy vitamin D<sub>2</sub>/D<sub>3</sub> were from 2.4 to 150 ng/mL. Typical calibration curves for analytes post extraction from stripped serum are shown in Fig. 2. The calibration curves that were generated had linear regression values of  $r^2$  >0.998 for each curve.



Fig. 2 Typical Calibration Curves for 25-OH Vitamin D<sub>2</sub>/D<sub>3</sub>

## Table 2 Comparison of ISOLUTE® SLE+ Treatment with Protein Precipitaion

25-hydroxy vitamin D2					ISOLUTE® SLE+ treatment					
Level	STD		Protein precipitation		Pre- spiked		Post- spiked		Matrix	Recovery
(ng/mL)	area	%RSD	area	%RSD	area	%RSD	Area	%RSD	Effect	
2.4	27,168	5.49	1,619	24.34	4,846	9.29	9,612	3.16	35 %	50 %
4.7	52,991	2.32	2,585	17.35	9,837	7.77	10,786	4.09	20 %	91 %
9.4	100,374	3.52	6,997	4.68	24,595	4.85	24,699	3.45	25 %	100 %
18.8	198,154	0.59	12,484	4.05	45,006	9.11	50,536	1.06	26 %	89 %
37.5	368,606	1.91	28,593	8.64	89,217	2.47	107,515	2.40	29 %	83 %
75	831,880	2.15	65,483	2.12	168,910	4.74	209,480	2.52	25 %	81 %
150	1,694,370	1.47	127,188	2.50	338,857	7.88	441,717	4.31	26 %	77 %
								Average	27 %	82 %

25-hydroxy vitamin D <sub>3</sub>					ISOLUTE <sup>®</sup> SLE+ treatment					
Level	STD		Protein precipitation		Pre- spiked		Post- spiked		Matrix	Recovery
(ng/mL)	area	%RSD	area	%RSD	area	%RSD	area	%RSD	Effect	
2.4	29,924	6.26	8,398	2.63	17,122	3.80	28,245	1.33	94 %	61 %
4.7	53,757	2.43	7,263	9.05	25,982	1.31	45,086	3.44	84 %	58 %
9.4	111,778	5.32	15,255	3.53	76,775	0.48	85,577	1.10	77 %	90 %
18.8	217,082	3.13	15,189	7.29	159,245	1.52	172,609	1.35	80 %	92 %
37.5	409,324	2.02	31,658	6.81	340,455	1.23	349,433	1.64	85 %	97 %
75	930,622	2.53	66,674	7.15	619,011	0.15	679,384	3.06	73 %	91 %
150	1,906,127	2.31	114,775	5.84	1,197,905	0.85	1,356,054	0.50	71 %	88 %
								Average	81 %	82 %

Recovery of analytes was determined by compare prespiked with post-spiked. The recoveries for analytes ranged above 80 % with RSDs <10 % (N=3).

Fig. 3 shows MRM chromatograms of 25-hydroxy vitamin  $D_2/D_3$  (each 4.7 ng/mL) treated by protein precipitation (A and B) or ISOLUTE<sup>®</sup> SLE+ (C and D). It is clear that ISOLUTE<sup>®</sup> SLE+ treatment provide a significantly more efficient extract than protein precipitation, leading to more reliable quantitation due to minimization of matrix effects.



Fig. 3 MRM Chromatograms of 25-OH Vitamin D<sub>2</sub>/D<sub>3</sub> (Each 4.7 ng/mL) Comparing ISOLUTE<sup>®</sup> SLE+ Treatment with Protein Precipitation

[Acknowledgement] We appreciate cooperation from Biotage Japan, Ltd..



For Research Use Only. Not for use in diagnostic procedures.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

www.shimadzu.com/an/

Shimadzu Corporation

First Edition: May. 2015