

# Application News



**Gas Chromatography** 

## Analysis of Residual Solvents in Pharmaceuticals by Water-Soluble Samples Using N<sub>2</sub> Carrier (JP17 Supplement II, USP 467)

The Japanese Pharmacopoeia 17<sup>th</sup> Edition (JP17) and the United States Pharmacopeia (USP) General Chapters <467> Residual Solvents provide test methods for residual solvents in pharmaceuticals mainly using headspace gas chromatography (GC). Residual solvents in pharmaceuticals are classified from Class 1 to 3 based on their potential human health risk. Since these compounds are strictly controlled, highly sensitive analysis is required. Helium (He) is generally used as the carrier gas, but as He supply shortages have become an issue recently, analysis using an alternative carrier gas such as N<sub>2</sub> has also been in demand lately.

This article introduces a JP17 Supplement II- compliant analysis of water-soluble samples for Class 1 and 2 residual solvents using N2 as the carrier gas.

N. Iwasa, T. Wada

## Instrument Configuration and Analysis Conditions

Class 1 and 2 standard solutions were prepared and measured in accordance with Procedure A and B of JP17 Supplement II, using Shimadzu HS-20 headspace gas sampler connected to Nexis<sup>™</sup> GC-2030 gas chromatograph. The two procedures differ in the type of column, the column temperature and the split ratio. Table 1 lists the GC and HS-20 analysis conditions used in this experiment.

GC analysis conditions (	,				
Model	: Nexis™GC-2030				
Detector	: FID-2030 flame ionization detector				
Column	: A) SH-Rxi™-624 Sil MS (0.32 mm l.D. × 30 m, d.f.= 1.8 μm)				
	B) SH-Stabilwax™ (0.32 mm l.D. × 30 m, d.f.= 0.25 μm)				
Column temp.	: A) 40 °C (20 min) - 10 °C /min - 240 °C (20 min) Total 60 mins				
	B) 50 °C (20 min) - 6 °C /min - 165 °C (20 min) Total 59.17 mins				
Injection mode	: A) Split 1:5				
	B) Split 1:10				
Carrier gas controller	: Linear velocity (N <sub>2</sub> )				
Linear velocity	: 35 cm/sec				
Detector temp.	: 250 °C				
FID H₂ flow rate	: 32 mL/min				
FID make up flow rate	: 24 mL/min (N <sub>2</sub> )				
FID air flow rate	: 200 mL/min				
Injection volume	: 1 mL				
HS-20 analysis conditions (same for Procedure A and B)					
Oven temp.	: 80 °C				
Sample line temp.	: 110 ℃				
Transfer line temp.	: 120 °C				
Vial stirring	: Off				
Vial volume	: 20 mL				
Vial heat-retention time	: 60 min				
Vial pressurization time	: 1 min				
Vial pressure	: 75 kPa				
Loading time	: 0.5 min				
Needle flush time	: 5 min				

## Analysis of Class 1 Standard Solution (Water-Soluble Sample)

The analysis results for Procedure A and B are shown in Fig. 1 and 2 respectively. Table 2 and 3 summarize the S/N ratios and the repeatability of Procedure A and B respectively.

The results obtained with Procedure A satisfied the requirements of JP17 Supplement II which states that "the S/N ratio for 1,1,1-trichloroethane in the Class 1 standard solution is not less than 5 and the relative standard deviation of each peak area is not more than 15 %." Satisfactory results were also obtained with Procedure B, as "the S/N ratio for benzene in the Class 1 standard solution is not less than 5 and the relative standard deviation of each peak area is not more than 15 %."

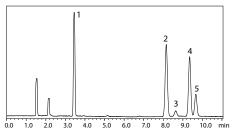


Fig. 1 Chromatogram of Class 1 Standard Solution by Procedure A (Water-Soluble Sample)

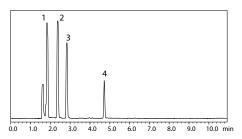


Fig. 2 Chromatogram of Class 1 Standard Solution by Procedure B (Water-Soluble Sample)

Table 2 S/N Ratio and Repeatability of Class 1 Standard Solution (Procedure A)

	•	•	
Peak No.	Compound	S/N ratio <sup>*1</sup>	%RSD (n=6)*1
1	1,1-Dichloroethane	263	1.81
2	1,1,1-Trichloroethane	242	1.74
3	Carbon tetrachloride	18	2.72
4	Benzene	204	2.08
5	1,2-Dichloroethane	69	1.56

#### Table 3 S/N Ratio and Repeatability of Class 1 Standard Solution (Procedure B)

Peak No.	Compound	S/N ratio <sup>*1</sup>	%RSD (n=6)*1
1	1,1-Dichloroethane	174	1.77
2	1,1,1-Trichloroethane	181	2.18
	+Carbon tetrachloride		
3	Benzene	142	1.12
4	1,2-Dichloroethane	105	1.02

In Table 2 and 3, the items specified in JP17 Supplement II are shown in red. \*1 The S/N ratios and relative standard deviation (%RSD) are reference values and not intended to be guaranteed values.

## Analysis of Class 2 Standard Solution (Water-Soluble Sample)


Fig. 3 and 4 below show the analysis results for Procedure A and B (Black: Class 2 mixture A standard solution (Class 2A), Pink: Class 2 mixture B standard solution (Class 2B), Blue: MIBK).

For system suitability, JP17 Supplement II specifies that "the resolution between acetonitrile and methylene chloride in the Class 2 mixture A standard solution is not less than 1.0" when using Procedure A, and "the resolution between acetonitrile and cis-1,2-dichloroethene in the Class 2 mixture A standard solution is not less than 1.0" when using Procedure B. Satisfactory results were obtained with both procedures.

\* The resolutions shown in the Fig. 3 and 4 are reference values and not guaranteed.

### Conclusion

In the analysis of residual solvents in water-soluble pharmaceuticals using the headspace GC method, satisfactory results comparable to those with He carrier gas were obtained with  $N_2$  as the carrier gas.

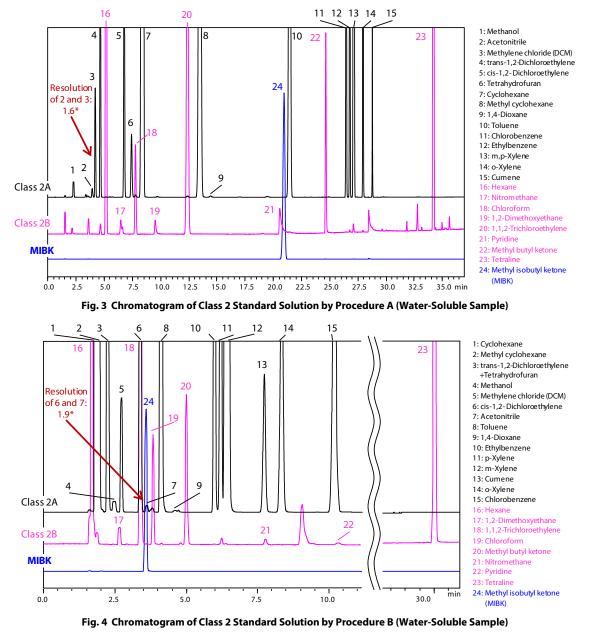

Nexis is a trademark of Shimadzu Corporation in Japan and/or other countries.

Rxi and Stabilwax are either trademarks or registered trademarks of Restek Corporation in the United States and/or other countries.

First Edition: Jun. 2020



Shimadzu Corporation www.shimadzu.com/an/

#### For Research Use Only. Not for use in diagnostic procedure.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. Shimadzu disclaims any proprietary interest in trademarks and trade names used in this publication other than its own. See <a href="http://www.shimadzu.com/about/trademarks/index.html">http://www.shimadzu.com/about/trademarks/index.html</a> for details.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.