

Application News

GC-MS GCMS-QP<sup>™</sup>2020 NX

# Fast Analysis of Fatty Acids in Brans by GC/MS

## No. M306


Y. Nakagawa

#### **User Benefits**

- Increased productivity with 10 mins run time for FAME 37
- EI/CI mode in the same batch to never misidentify a target FAME peak
- Constant linear velocity mode to easily convert the method to GC-FID

#### Introduction

Fatty acids, particularly those with a high degree of unsaturation

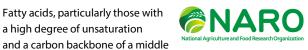

chain length, are known as functional nutrients.

Shimadzu Corporation Japan has entered the second year of the three year collaborative research with National Agriculture and Food Research Organization (NARO) on the amount of functional nutrients (e.g. fatty acids) in foods such as brans, tea leaves and rice.

As a part of the collaboration, fatty acids (Fig. 2) were quantitated in 48 bran samples. Details of the results are off limits to the public at the moment of this writing.

However, the otherwise time-consuming run of the 48 samples was cut to 10 mins by optimizing a method without using a high heat ramp rate, which warranted this application news.



Fig. 1 GCMS-QP<sup>™</sup>2020 NX

#### Methods - Extraction

Extraction was performed according to Fatty Acid Methylation Kit (Nakarai USA Inc., P/N MAK224-1KT) with the sample weight of 0.1 ±0.005 g.

### Methods – Analysis

Table 1 Instrument Configurations		
GC-MS	: GCMS-QP 2020 NX	
Auto Injector	: AOC™-20i Plus	
Auto Sampler	: AOC-20s Plus	
Analytical Column	: DB-FastFAME (20 m $\times$ 0.18 mm l.D., df=0.20 $\mu$ m) P/N: G3903-63010	
Glass Insert	: Split liner with wool	

Table 2 Analytical Conditions		
GC		
Inlet temp.	: 250 °C	
Injection Mode	: Split	
Spit ratio	: 10	
Carrier gas	: Helium	
Control Mode	: Constant linear velocity (53.4 cm/s)	
Column oven temp.	: 60 °C (1 min) $\rightarrow$ (40 °C /min) $\rightarrow$ 200 °C (3 mins) $\rightarrow$ (25 °C /min) $\rightarrow$ 250 °C (1 min) Total 10.50 mins	
Purge flow rate	: 5 mL/min	
Sample Inj. volume	: 1 μL	
MS		
Ion Source Temp.	: 230 °C	
Interface temp.	: 250 °C	
CI reagent gas	: Isobutane	
Measurement Mode	: Simultaneous Scan/SIM (FAAST)	
Scan mass range ( <i>m/z</i> )	: 35-600 at scan speed of 20000 $\mu/s$	

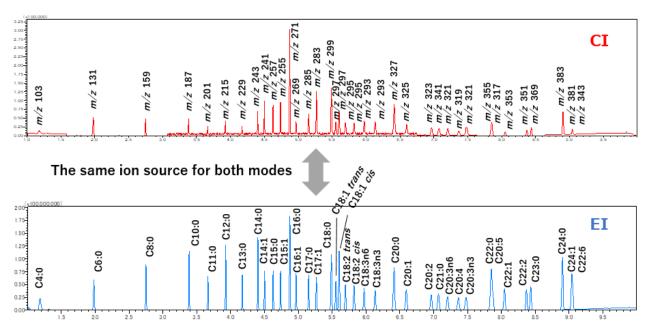


Fig. 2 CI and El Chromatogram at 20 µg/mL with Smart Ion Source

#### Smart EI/CI Ion Source




It is of paramount importance to have CI spectra for FAME analysis. CI spectra helps determine elution positions of FAMEs that might vary due to column lot difference and column conditions on the day of analysis.


Smart EI/CI ion source allows a seamless switching of EI and CI ionization modes within the same batch (Fig. 2). When the identity of a peak is questioned, the CI data from the same batch, as opposed to comments such as "retention time won't change" or "vendor catalog said this peak is this", will satisfy customers and auditors alike.

The *m/z* in Fig. 2 are quasi-molar masses for illustration and not necessarily those obtained.

#### Calibration Curve

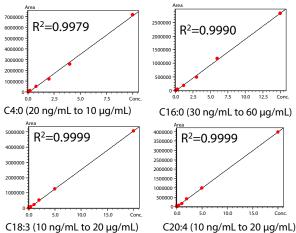

#### Importance of CI Mode

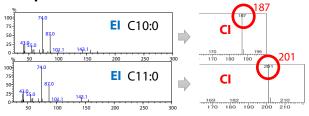

### Summary

Fast analysis of fatty acids in brans were performed with Smart El/Cl ion source. Cl data was acquired in the same batch without changing the ion source. It is of utmost importance to have CI data ready at hand to ensure the confidence in peak identity on the day of analysis, given the variability due to the column lot and column conditions.

Calibration curves for Fame 37 were all satisfactory with R<sup>2</sup> of greater than 0.990. 48 bran samples, the results of which are in non-disclosure agreement, were analyzed with the calibration curves

Shimadzu Corporation would like to thank Hironori Juichi, Yayoi Ichiki and Mari Yamamoto from NARO for their guidance in leading the project.

GCMS, GCMS-QP, GCMS-TQ and AOC are trademarks of Shimadzu Corporation in Japan and/or other countries.



#### Shimadzu Corporation

Analytical & Measuring Instruments Division **Global Application Development Center** 

www.shimadzu.com/an/

First Edition: Dec. 2020

For Research Use Only. Not for use in diagnostic procedure. This publication may contain references to products that are not available in your country. Please contact us to check the availability of these

products in your country. The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. See http://www.shimadzu.com/about/trademarks/index.html for details. Third party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they

The information contained here in is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or radius

completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.