

Detection of Reductively-Alkylated Cysteines Using Protein Sequencer

Introduction

The main stream in current protein analysis is proteome analysis utilizing a mass spectrometer and a search engine based on genomic databases to identify proteins. Proteins form specific 3-dimensional structures and display the functions of the protein concerned by having that 3-dimensional structure. The 3-dimensional structures of proteins consist of four levels, one of which is the tertiary structure formed by pairs of the side chains of the amino acids that comprise the protein. One of these 3-dimensional structures is created by intramolecular crosslinking by disulfide (S-S) bonds of cysteines in a molecule. Disulfide bonds have the function of stabilizing the 3-dimensional structure of the protein and normally form S-S bonds in many cases. For this reason, it is sometimes necessary to identify cysteines when conducting an amino acid sequence analysis. When using a protein sequencer, even if the half-cystine residue on N-terminal side is converted to ATZ-half-cystine by Edman degradation, the half-cystine residue will form a disulfide bond with the half-cystine on C-terminal side, and elution and detection by HPLC will be impossible. Therefore, sequence analysis is generally carried out after reductive alkylation of the protein to sever the disulfide bonds, and the cysteine is then identified as a PTH-modified cysteine.

This article introduces an example of identification of a PTH-modified cysteine (phenylthiohydantoin derivative of reductively-alkylated cysteine) by analyzing reductively-alkylated peptides/proteins with PPSQ™ -50A series protein sequencer systems (isocratic system and gradient system).

T. Kuriki

Reductive Alkylation Protocol

Several reductive alkylation methods for samples containing disulfide bonds are available. Here, dithiothreitol (DTT) was used as the reductant, and alkylation was conducted by carboxymethylation using monoiodoacetic acid, carbamidomethylation using iodoacetamide, and pyridylethylation using 4-vinylpyridine. Fig. 1 shows the reaction protocol of reductive alkylation. In order to apply the treatment samples directly to the PPSQ-50A series without purification, N-ethylmorpholine was used in the reaction solution, as it has little effect on Edman degradation. Because the optimum pH differs depending on the alkylation agent, pH was adjusted by using an appropriate concentration of N-ethylmorpholine.

The sample used here was 100 pmol synthetic oxytocin containing one intramolecular disulfide bond (Peptide Institute, Inc., Code: 4084-v).

Table 1 and Table 2 show the conditions of the analysis by the protein sequencer when using the isocratic system and the gradient system, respectively.

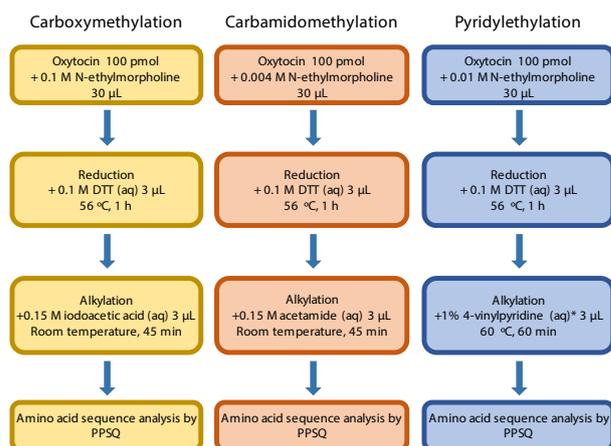


Fig. 1 Reaction Protocol of Reductive Alkylation

Table 1 Analysis Conditions (Isocratic System)

Column	: Wakopak™ Wakosil® PTH-II (250 mm × 4.6 mm I.D.)
Mobile phase	: PTH-amino Acids Mobile Phase
Flow rate of mobile phase	: 1.0 mL/min
Column temp.	: 40 °C
Detection	: SPD-M30A (269 nm) with High Sensitivity Flow cell

Table 2 Analysis Conditions (Gradient System)

Column	: Wakopak™ Wakosil® PTH-GR (S-PSQ) (250 mm × 2.0 mm I.D.)
Mobile phase	: A: PTH-amino Acids Mobile Phase A (for Gradient Elution) B: PTH-amino Acids Mobile Phase B (for Gradient Elution)
Flow rate of mobile phase	: 0.3 mL/min
Column temp.	: 35 °C
Detection	: SPD-M30A (269 nm) with High Sensitivity Flow cell

Next, 3 µL (30 pmol) of the reductively-alkylated sample solution was added to a PVDF membrane and dried. After drying, a further polybrene solution was added to the PVDF membrane and the sequence analysis was started. Fig. 2 and Fig. 3 show the sequence analysis results for oxytocin. Fig. 3 is an enlarged view of part of Fig. 2.

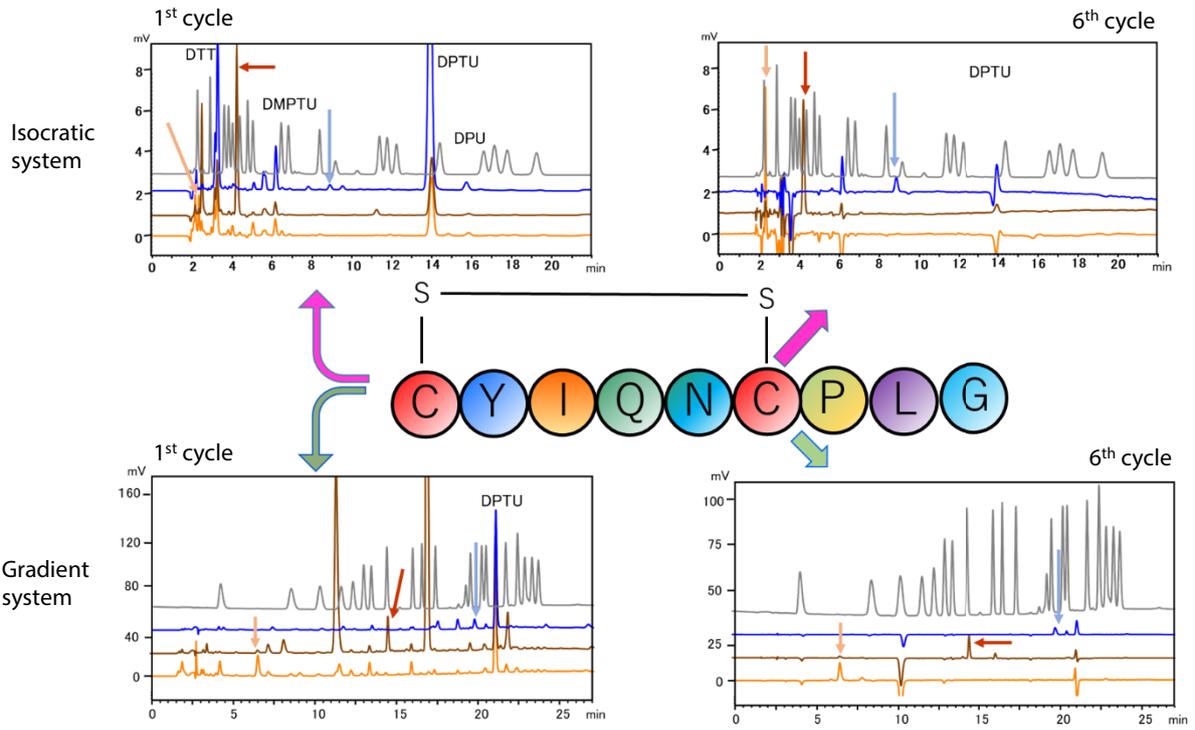


Fig. 2 Amino Acid Sequence of Reductively-Alkylated Oxytocin and Its Chromatograms
Orange: Carboxymethylation (CM), Brown: Carbamidomethylation (CAM), Blue: Pyridylethylation (PE);
1st Cycle : Raw Chromatogram, 6th Cycle : Subtraction Chromatogram

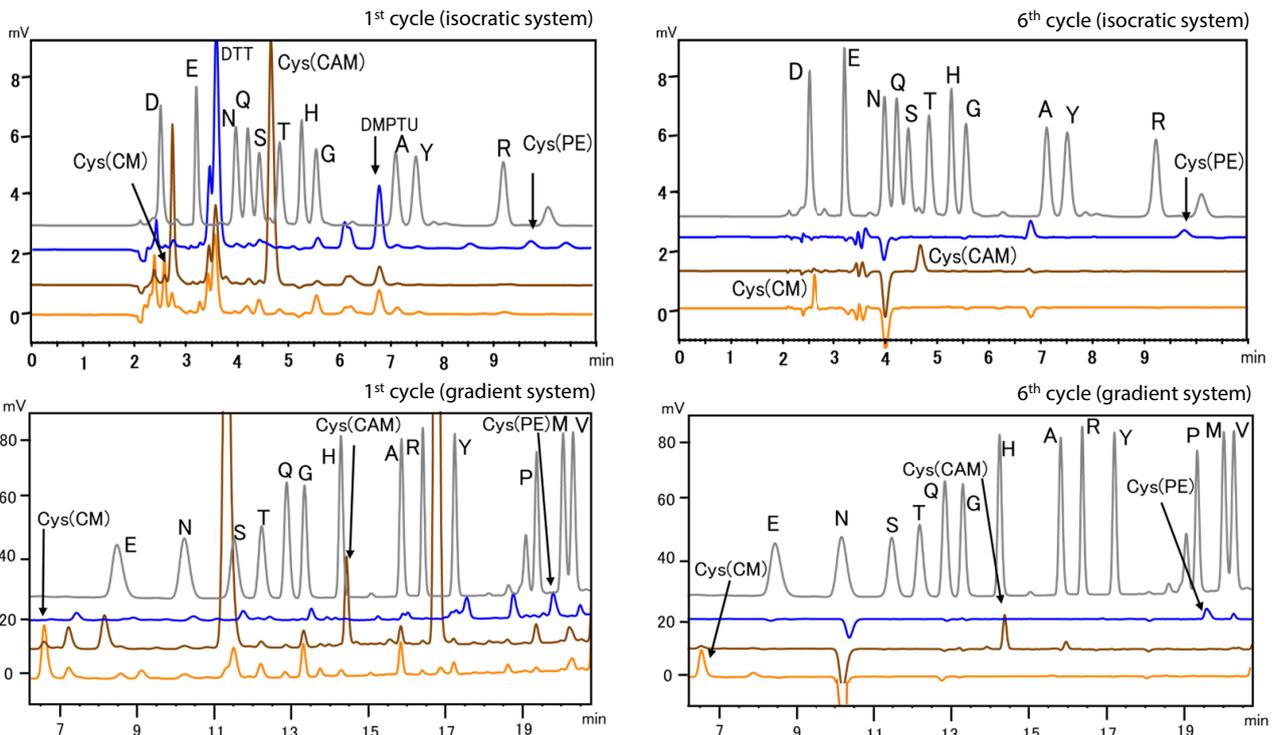


Fig. 3 Amino Acid Sequence of Reductively-Alkylated Oxytocin and Its Chromatograms (Partial Enlargement of Fig. 2)
Orange: Carboxymethylation (CM), Brown: Carbamidomethylation (CAM), Blue: Pyridylethylation (PE);
1st Cycle : Raw Chromatogram, 6th Cycle : Subtraction Chromatogram

From these analysis results, when using the isocratic system, elution of the cysteine occurred in the vicinity of 2.4 min with PTH-carboxymethylation, 4.3 min with PTH-carbamidomethylation, and 8.9 min with PTH-pyridylethylation. However, because the cysteine was detected near the elution position of PTH-Asp with PTH-carboxymethylation and near that of PTH-Thr with PTH-carbamidomethylation, care is necessary in identification. In contrast, identification is easy with pyridylethylation, as elution of the PTH-pyridylethylcysteine occurred at 8.9 min and did not overlap with the other PTH-standard amino acids. On the other hand, when using the gradient system, elution of the cysteine occurred in the vicinity of 6.5 min with PTH-carboxymethylation, 14.3 min with PTH-carbamidomethylation, and 19.7 min with PTH-pyridylethylation. With PTH-carbamidomethylcysteine, detection occurred near the elution position of PTH-His, indicating that care is required in identification. Essentially, preparation of protein sequencer samples by desalination is recommended in order to avoid salt contamination of the samples as far as possible. In this report, it was possible to conduct the sequence analyses without desalination by reducing the concentrations of salts that affect Edman degradation. However, in sample preparation by reductive carbamidomethylation, peaks originating from reductive alkylation reagents were detected in the vicinity of 11.2 min and 16.6 min (Figs. 2 and 3). Desalination treatment of the samples is necessary in order to remove these peaks.

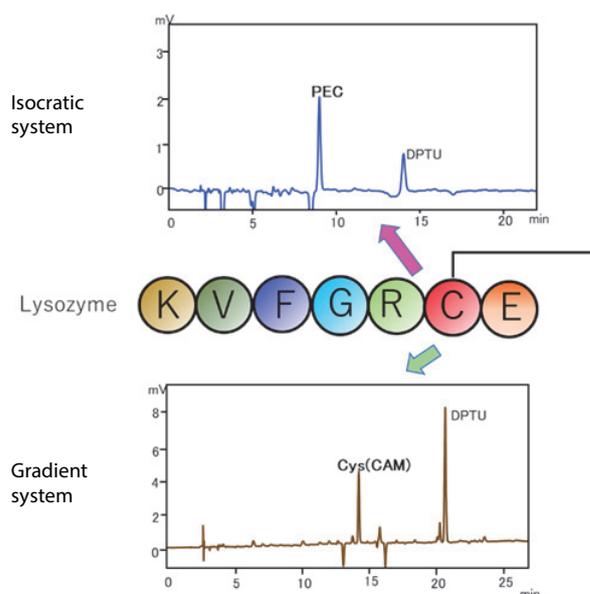


Fig. 5 6th Cycle Chromatograms of Lysozyme after Reductive Alkylation and Edman Degradation in Reactor (Subtraction Chromatograms)
PEC: Pyridylethylcysteine
Cys(CAM): Carbamidomethylcysteine

Reductive Alkylation of Cysteine Residue in Protein Sequencer

To facilitate identification of cysteine residues, Shimadzu PPSQ-50A series protein sequencers include an analysis program that conducts reductive alkylation in the instrument, followed directly by Edman degradation. The isocratic system uses a reaction program in which pyridylethylation of the sample is conducted in the reactor, while the gradient system uses a reaction program for sample carbamidomethylation in the reactor. Fig. 4 shows the respective reaction schemes. Fig. 5 shows the chromatograms for the 6th cycle of a sequence analysis of 30 pmol of lysozyme (Sigma-Aldrich Co. LLC, product No. L6876). Cysteine residue which has formed a disulfide bond with C-terminal cysteine residue exists in the 6th cycle, demonstrating that cysteine residues which have been stably modified by the dedicated reaction program can be detected easily.

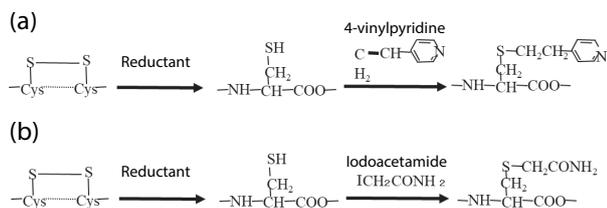


Fig. 4 Reaction Schemes
(a) Pyridylethylation, (b) Carbamidomethylation

PPSQ is a trademark of Shimadzu Corporation in Japan and/or other countries.
Wakopak and Wakosil are either registered trademarks or trademarks of FUJIFILM Wako Pure Chemical Corporation.

Conclusion

In comparison with other amino acids, sample preparation by reductive alkylation to be detected stably is necessary in analyses of cysteines and cystines, which decompose during Edman degradation if analyzed without appropriate pretreatment, and a certain amount of the sample is also necessary for identification.

The Shimadzu PPSQ-50A series protein sequencers enable easy and accurate identification of N-terminal sequence. In addition, easy identification of unmodified cysteines and cystines is possible, as reported in Application News No. B111, and identification of the reductively-alkylated cysteine residues in this report is also easy. Thus, the PPSQ-50A protein sequencers provide an effective tool for structural analysis in proteome analysis and for evaluation of synthetic intermediates before the formation of disulfide bonds in peptide synthesis.

First Edition: Nov. 2020



Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedure.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. Shimadzu disclaims any proprietary interest in trademarks and trade names used in this publication other than its own. See <http://www.shimadzu.com/about/trademarks/index.html> for details.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

© Shimadzu Corporation, 2020