

Application News

No.042

Total Organic Carbon Analysis

Recovery Test for Sodium Dodecylbenzenesulfonate by TOC-L_{CPH}

In the test method for total organic carbon as listed in the General Tests, Processes and Apparatus section of the Japanese Pharmacopeia (JP), it is specified that the TOC analyzer must be capable of measurement of sample concentrations below 0.050 mg/L, and that the instrument achieve a recovery rate of at least 0.450 mg/L carbon in measurement of a sodium dodecylbenzenesulfonate solution (0.806 mg/L). Here, we introduce an example of recovery testing with a sodium dodecylbenzenesulfonate solution using the Shimadzu combustion catalytic oxidation model TOC-LCPH, and an example of analysis of a sample with a TOC concentration below 0.050 mg/L.

Recovery Testing for Sodium Dodecylbenzenesulfonate

The Japanese Pharmacopeia (JP) specifies that "the decomposition device for either method should be capable of generating not less than 0.450 mg/L of organic carbon when using a solution of sodium dodecylbenzenesulfonate (theoretical value of total organic carbon in this solution is 0.806 mg/L) as the sample."

Because sodium dodecylbenzenesulfonate solution (0.806 mg/L) has a carbon concentration of 0.500 mg/L, the instrument must be capable of combustion oxidation sufficient to permit a 90 % detection rate (recovery) when measuring this solution.

Thus, we prepared a sodium dodecylbenzenesulfonate solution with a carbon concentration of 0.500 mg/L by dissolving sodium dodecylbenzenesulfonate in distilled water, and then conducted TOC measurement (TC-IC measurement). The results are shown in Fig. 1 and Table 1.

We calibrated the instrument by generating calibration curves using 0 and 1 mgC/L (1 mg/L carbon concentration) standard solutions for both the TC and IC measurements. To eliminate the influence of carbon in the distilled water used to prepare the standard solutions for generating both calibration curves, we performed correction using the calibration curve Shift to Origin feature.

Measurement of the sodium dodecylbenzenesulfonate solution with a carbon concentration of 0.500 mg/L yielded a result of 0.498 mg/L (498 µg/L), which is a recovery rate of 99.6 %, thus satisfying the instrument specifications stipulated in the total organic carbon test method of the Japanese Pharmacopeia (JP).

<measurement con<="" th=""><th>ditions></th></measurement>	ditions>
Instrument	: Shimadzu TOC-LCPH total organic
	carbon analyzer
Measurement Item	: TOC (TOC by TC-IC)
Calibration Curves	: TC; 0 to 1 mgC/L using potassium
	hydrogen phthalate aqueous
	solutions
	IC; 0 to 1 mgC/L using sodium
	carbonate/sodium bicarbonate
	aqueous solutions
Sample	: Sodium dodecylbenzenesulfonate
1	(Wako Pure Chemical Industries,
	Ltd.) reagent (for ABS
	measurement) was dissolved in
	distilled water to prepare a
	0.806 ma/L solution (0.500 ma/L
	carbon concentration).



Fig. 1 TOC Measurement Data for 0.500 mg/L (500 µg/L) Sodium Dodecylbenzenesulfonate Solution

Table 1 TOC Measurement Data for 0.500 mg/L (500 µg/L) Sodium Dodecylbenzenesulfonate Solution

Sample Name	TC Value [mgC/L]	IC Value [mgC/L]	TOC (TC-IC) Value [mgC/L]	Recovery [%]
Sodium dodecylbenzenesulfonate solution with a carbon concentration of 0.500 mg/L	0.552	0.054	0.498	99.6

TOC Measurement Below 0.050 mg/L

The Japanese Pharmacopeia (JP) specifies that "the instrument should be capable of measuring the amount of organic carbon down to 0.050 mg/L."

Thus, we prepared a potassium hydrogen phthalate aqueous solution with a TOC concentration of 0.025 mgC/L. The results are shown in Fig. 2 and Table 2. Because the distilled water used to prepare the sample includes some slight amount of organic carbon-containing substances, the result of measurement of the 0.025 mgC/L potassium hydrogen phthalate aqueous solution was 0.0407 mgC/L, but the coefficient of variation (CV value) was 2.42 %. This result is near the lower limit of quantitation, and within 10 % of the CV value, demonstrating that the Shimadzu TOC-LCPH combustion catalytic oxidation type TOC analyzer satisfies the TOC measurement criteria value of less than 0.050 mg/L stipulated in the Japanese Pharmacopeia total organic carbon test method.



Fig. 2 Data Obtained with Less than 0.05 mg/L (50 µg/L) TOC

<measurement conditions=""></measurement>				
Instrument	:Shimadzu ТОС-L _{СРН}			
	combustion catalytic oxidation			
	type TOC analyzer			
Measurement Method	: TOC (TOC measurement using			
	acidification and sparging method)			
Calibration Curve	: 0 to 0.5 mgC/L using potassium hydrogen phthalate			
Injection Volume				
Sample	: 0.025 mgC/L notassium			
Sample	hydrogen phthalate aqueous solution			

Table 2 Measurement Data Obtained with Less than 0.05 mg/L (50 µg/L) TOC

Sample Name	TC Value [mgC/L]	CV Value [%]	
Potassium hydrogen phthalate aqueous solution	0.0407	2.42	

First Edition: February, 2012



For Research Use Only. Not for use in diagnostic procedures.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

Shimadzu Corporation www.shimadzu.com/an/