

Affinity purification of IdeZ digest for Glycosylation profile of Immunoglobulins using a linear benchtop MALDI-TOFMS.

ASMS 2019 WP-053

Y. Yamazaki¹, S. Nakaya¹
1 Shimadzu Corporation, Kyoto, Japan

Affinity purification of IdeZ digest for Glycosylation profile of Immunoglobulins using a linear benchtop MALDI-TOFMS.

Overview

- A batch analysis of glycosylation of IgG using a bench-top linear MALDI-TOFMS was intended.
- A magnet beads linked with protein A was applied to affinity purification, following a digestion of IgG by IdeZ.
- Three different glycosylated Fc were clearly observed to classify them with a statistical software.

Introduction

Glycosylation on protein plays wide-range vital roles in biological processes from stabilization of protein conformation to expression of binding specificity. In this view, a characterization of the N-/O-linked glycan is quite significant, especially, in development of biopharmaceuticals. To date, whereas intensive efforts were conducted to characterize glycans precisely with high-end mass spectrometers, conventional instruments without time consuming preparation has been anticipated

for batch analysis in screening or QA/QC. A newly developed bench-top MALDI-TOFMS is expected to be the conventional instrument in terms of sufficient specification, through-put, and cost effectiveness. We attempted to characterize glycosylation of IgG without a release of glycan using the bench-top MALDI-TOFMS. To do this, we examined a preparation using affinity beads and enzymatic cleavage.

Methods

• IgGs and their preparations

NIST mAb, IgG from mouse serum, and myeloma IgG were subject to MS analysis. IdeZ and Protein-A magnetic beads (Mag Sepharose) were purchased from Promega, and GE Healthcare respectively. The IgGs dissolved in tris-NaCl buffer were incubated with IdeZ at 37°C for two hours. Resultant solutions including Fc and F(ab')₂ were applied to the beads equilibrated with the tris-buffer, then washed with the same buffer several times. Fc associated with Protein-A was eluted with acidic solution, then desalted with ZipTip C18.

• MALDI-TOFMS

All MS analysis in positive ion mode was performed with a bench-top linear MALDI-TOFMS (MALDI-8020, Shimadzu Corp., Japan). Sinapinic acid (SA) and Ferulic acid (FA) were applied to MS analysis as matrix at the same condition as follows; 20 mg/mL in 50% acetonitrile/milliQ water with 0.1% TFA.

• Statistical data analysis

eMSTAT Solution (Shimadzu Corp., Japan) was applied to perform classification/differentiation of glycosylated-Fc regions.

Affinity purification of IdeZ digest for Glycosylation profile of Immunoglobulins using a linear benchtop MALDI-TOFMS.



MALDI-8020

Features:

- Compact design
- > 5,000 MS resolution
- High through-put with solid state laser and fast stage motion
- Rapid sample introduction
- UV laser-based source cleaning
- Quiet operation

Results

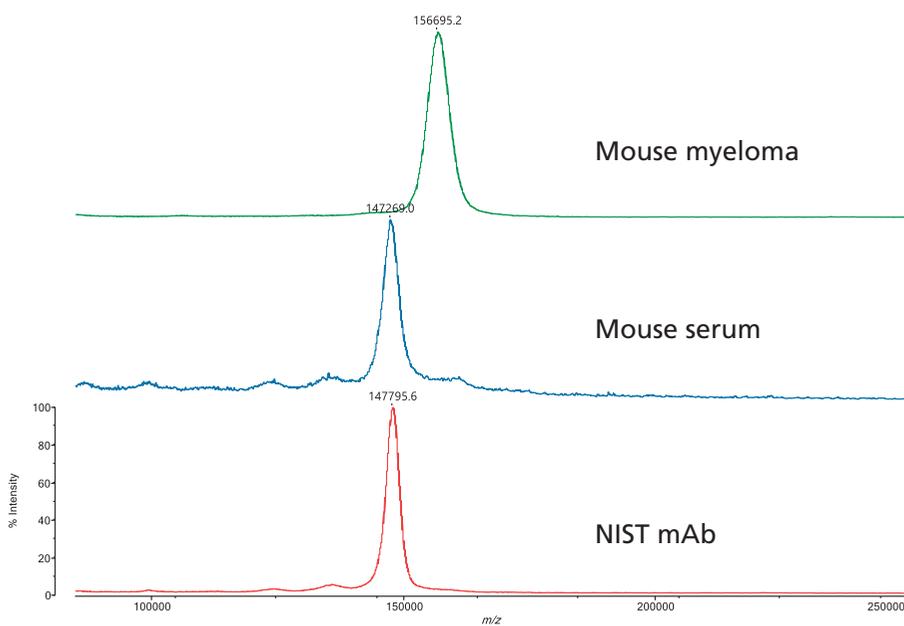


Fig.1 MS of whole IgGs

Affinity purification of IdeZ digest for Glycosylation profile of Immunoglobulins using a linear benchtop MALDI-TOFMS.

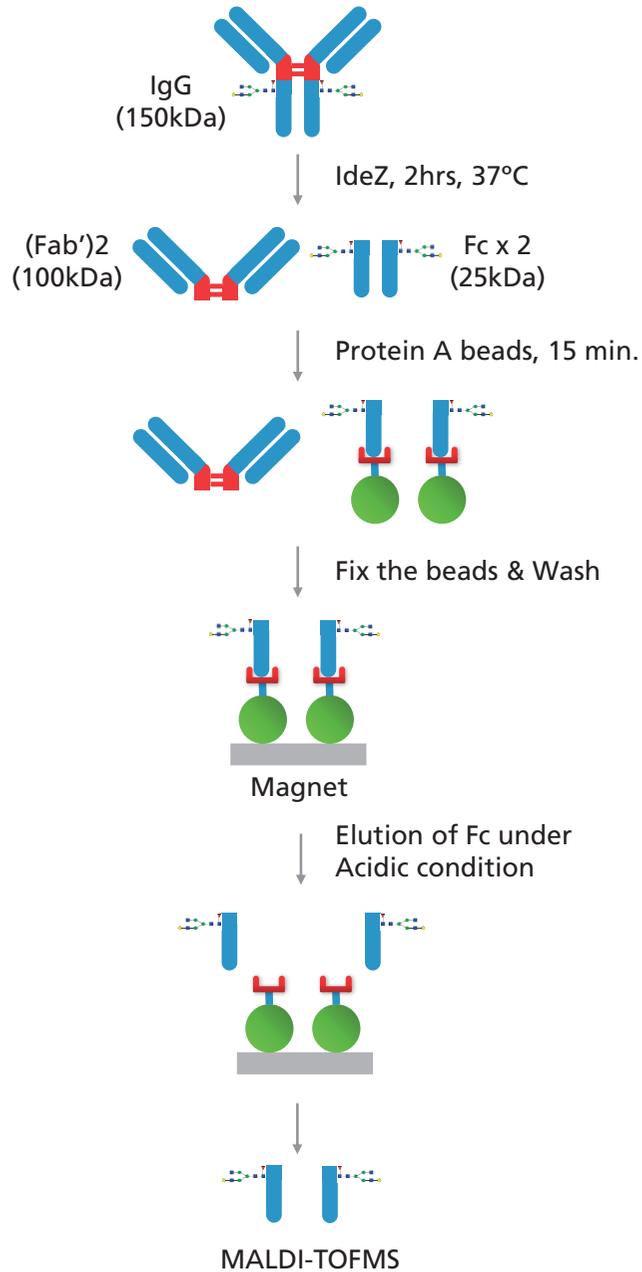


Fig.2 Work-flow of Fc affinity separation/detection

Affinity purification of IdeZ digest for Glycosylation profile of Immunoglobulins using a linear benchtop MALDI-TOFMS.

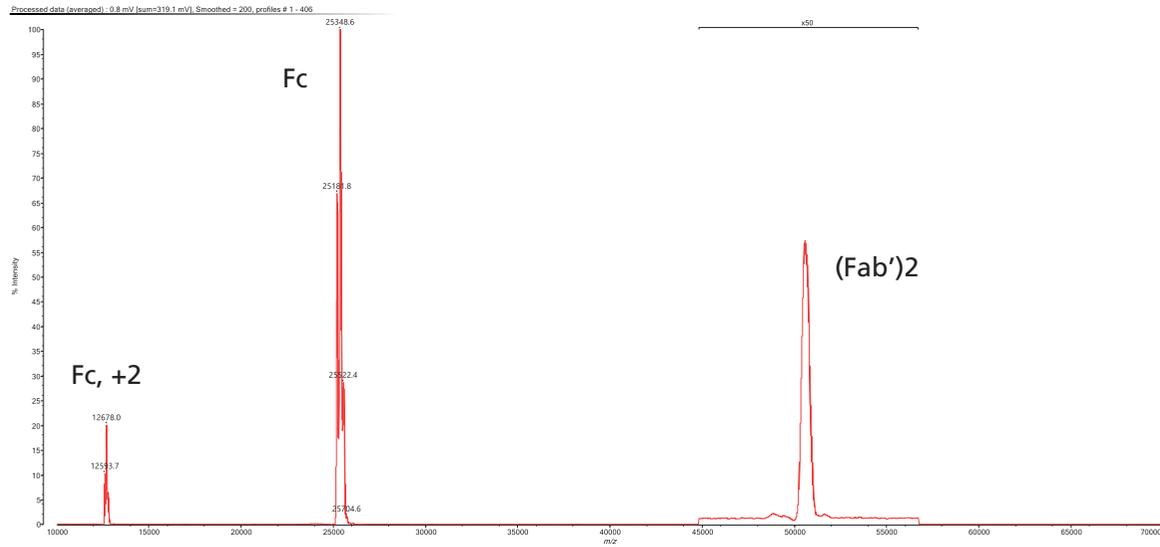


Fig.3 MS of NIST mAb treated with IdeZ

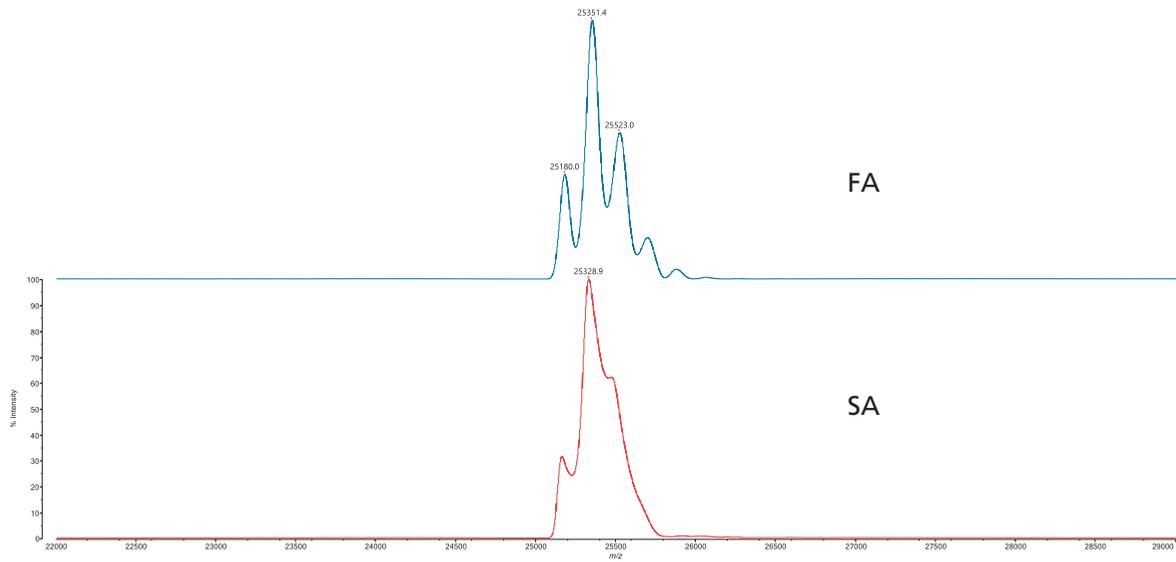


Fig.4 A comparison of Fc measured with SA and FA under the identical MS condition

Affinity purification of IdeZ digest for Glycosylation profile of Immunoglobulins using a linear benchtop MALDI-TOFMS.

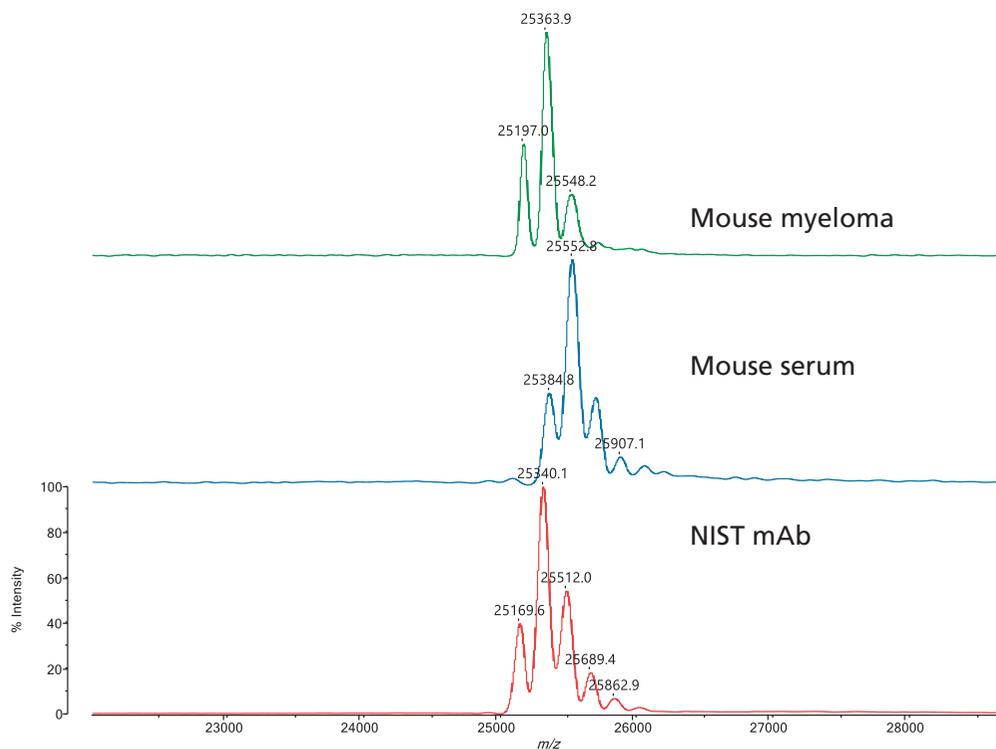


Fig.5 A comparison of three Fc separated with the work-flow

Affinity purification of IdeZ digest for Glycosylation profile of Immunoglobulins using a linear benchtop MALDI-TOFMS.



Fig.6 Classification of three Fc using eMSTAT Solution

Affinity purification of IdeZ digest for Glycosylation profile of Immunoglobulins using a linear benchtop MALDI-TOFMS.

Conclusions

- Purification of enzymatically fragmented IgGs by affinity-beads enables batch analysis for the glycosylation using a bench-top MALDI-TOFMS.
- MS resolution of the bench-top MALDI-TOFMS is sufficient to recognize three Fc that differ in glycosylation mainly.
- A statistical analysis by eMSTAT Solution enables a classification of three glycosylated Fc smoothly and quickly, which could be applicable to QA/QC.

For Research Use Only. Not for use in diagnostic procedures. This presentation may contain references to products that are not available in your country. All rights reserved. Information subject to change without notice.

First Edition: May, 2019