

Improving Preparative Scale Fractionation


Fraction Collector FRC-40





C193-E033A

Efficient, Easy to Use, Space Saving

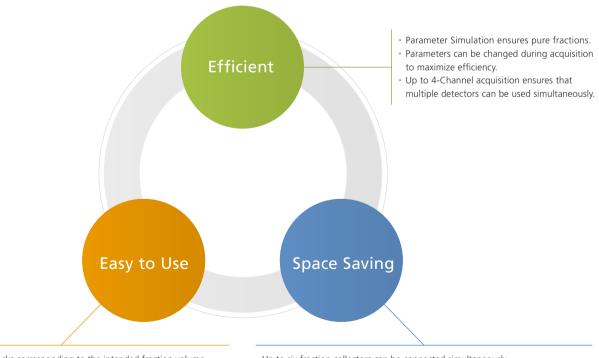

Compatible with many racks corresponding to the intended fraction volume.

Open access design allows a single system to be shared by multiple users.
Innovative sample rescue function minimizes sample loss when errors occur.

• Up to six fraction collectors can be connected simultaneously.


· Can accommodate up to 3,240 fractions (6 FRC-40s, 10 mm tubes).



Improved Purification Workflow

The FRC-40 incorporates greatly improved fraction simulation. Simply select the peaks of interest and the software dynamically configures parameters that will collect them. (Patent pending)

System parameters such as the pump flow rate and detector settings can be changed during analysis. This allows the user to fractionate during separation development and reduces the potential of wasting precious sample.

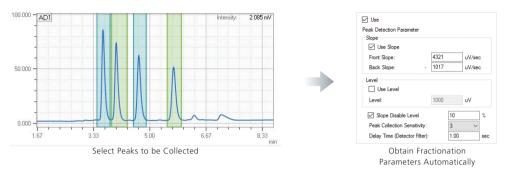

Simplify Purification Procedures

The FRC-40 is compatible with a variety of MTPs, vials, test tubes and bottles. This allows users to select the fraction size that best fits their analysis and maximizes the number of fractions that can be collected.

Racks are specified through the software, allowing multiple users to collect fractions without the risk of cross-contamination or mix-up. The open access design of the FRC-40 allows racks to be removed and replaced during analysis. This allows multiple purifications to take place using a single HPLC system.



Standard Fraction Containers* 10, 18, 25, and 35mm test tubes 96-well MTP/ DWP 250, 500, and 1,000 mL bottles

*Some custom containers may be used. Inquire if interested.



Use Bench Space Effectively

The instrument footprint has been optimized to accommodate the maximum number of fractions in the minimum amount of space. Space savings of up to 50% can be achieved compared to competing products.

Up to six FRC-40 units may be multiplexed to a single HPLC system, allowing users to customize the instrument capacity to their needs.

	Maximum Capacity (1 x FRC-40) / (6 x FRC-40)
10 mm test tubes 18 mm test tubes 25 mm test tubes 35 mm test tubes 250 mL bottle 500/1,000 mL bottles	540 / 3,240 216 / 1,296 108 / 648 54 / 324 20 / 120 12 / 72
96-well MTP/ DWP	9/ 54



LabSolutions is a trademark of Shimadzu Corporation.



Shimadzu Corporation www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures. This publication may contain references to products that are not available in your country. Please contact us to check the availability of these

products in your country. Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®". Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "@". Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The contents of this publication are provided to you "as is" without warranty of any kind, and are subject to change without notice. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication.