IMAGEREVEAL MS

Install Manual

Read this manual thoroughly before you use the product.
Keep this manual for future reference.

1 Install Manual (ENG)

1.1 Installing and Uninstalling

1.1.1 Installing

This section explains the procedure for installing all IMAGEREVEAL programs to your computer.

Use the same procedure when upgrading the version of your IMAGEREVEAL to overwrite the installation.

NOTE Depending on your anti-virus software, the installation may fail. In this case, disable your anti-virus software when installing this software. When upgrading IMAGEREVEAL, be sure to exit the software before installing.

1

Install MATLAB Runtime.

- 1 Unzip the [MATLABRuntime] folder on the installation disk.
As a preliminary step, create an empty extraction destination folder anywhere on the PC you are installing.

NOTE MATLAB must be installed before installing IMAGEREVEAL.
If the correct version of MATLAB Runtime is already installed, this procedure is unnecessary.


- 2 Select the [MATLABRuntime] folder on the installation disk and select [Expand All...] from the right-click menu.
The destination folder selection screen appears.
- 3 Set the extraction destination folder that you created in advance for the destination.
- 4 Click [Extract].
The extraction process is performed.

NOTE The extraction process may take several minutes to complete.

- 5 When the extraction is complete, double-click "setup.exe" in the extraction destination folder.
Start the MATLAB installer.


NOTE It may take a few minutes for the installation window to appear.

- 6 Click [Next].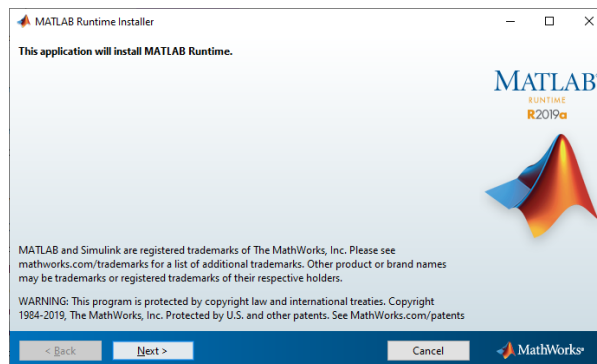


MATLAB Installation Window

- 7 Check the license agreement. If you agree, check the [Yes] checkbox and click [Next].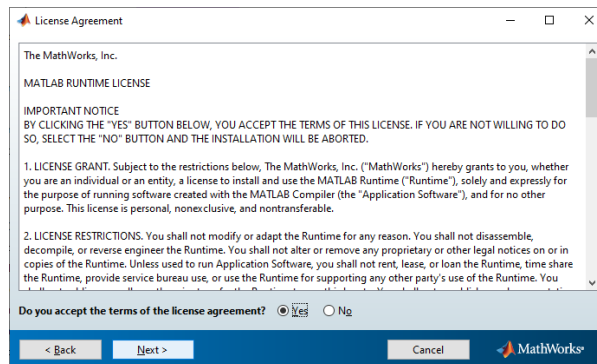

MATLAB License Agreement


8 Leave the installation folder as-is and click [Next].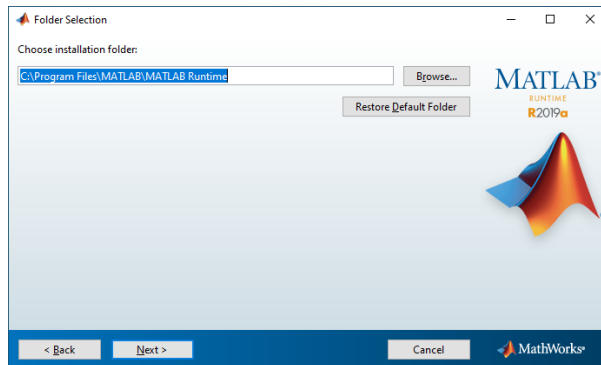

MATLAB Folder Selection

NOTE Do not change the installation folder from the default.

9 Confirm the settings and click [Install].

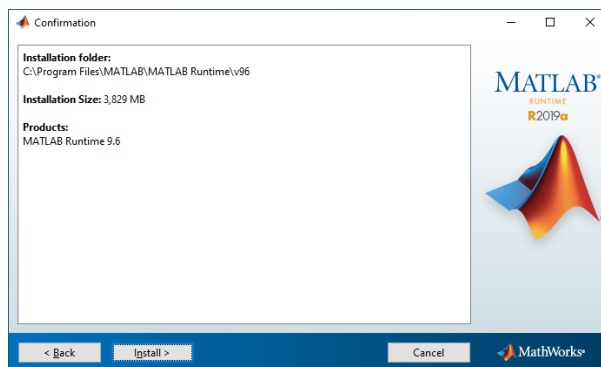
MATLAB Confirmation Window

The installation starts.

MATLAB Installation Progress Window

- 10 After the installation ends successfully, click [Exit] to close the window.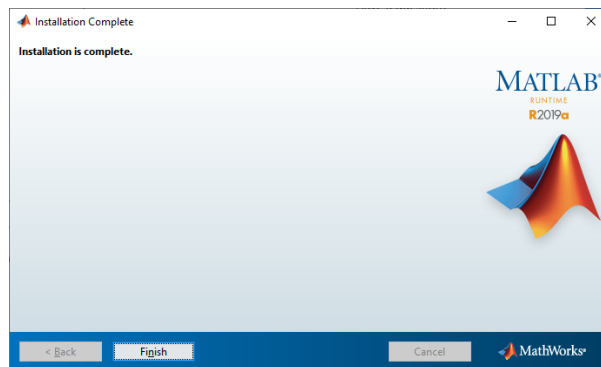

MATLAB Installation Complete Window


2

Install IMAGEREVEAL.

- 1 Double-click "IMAGEREVEALSetup.exe".


NOTE If MATLAB Runtime is not installed, IMAGEREVEAL cannot be installed.

- 2 Check the license agreement and conditions. If you agree, check the checkbox and click [Install].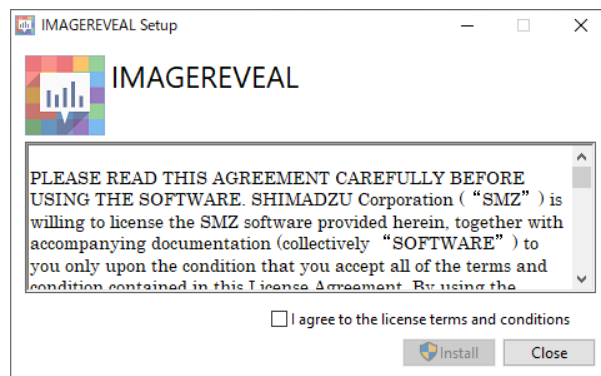


Installation Window

The installation starts.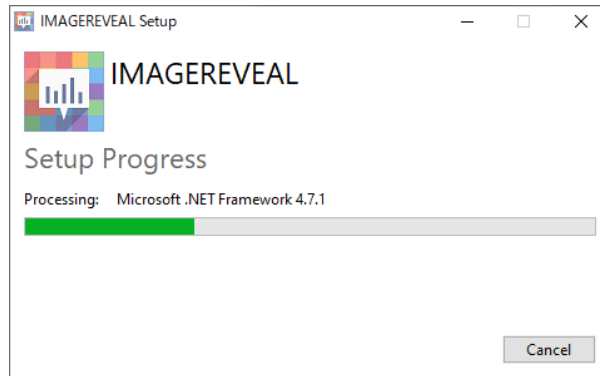

Installation Progress Window


NOTE If the following window is displayed, click [Yes].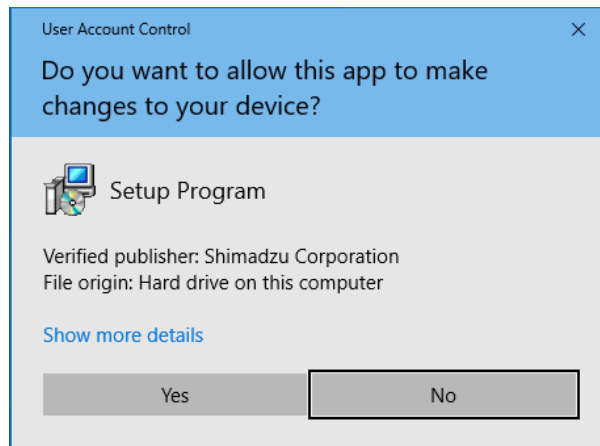

[User Account Control] Window

3

When the installation complete window is displayed, click [Restart] to restart your computer.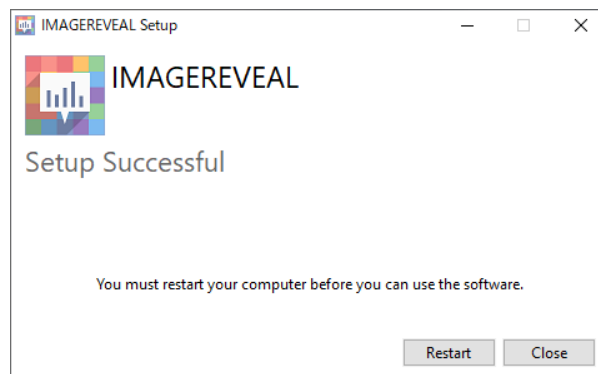

Installation Complete Window

NOTE

- If you have installed anti-virus software that performs real-time scans when files are accessed, it may take more time to perform operations such as saving the project file or creating a data matrix. Configure the anti-virus settings so that the folders and files related to this software are exempt from the scans.
- After this software is installed for the first time, you can use it as a 30-day demo version. However, the demo version has the following functional restrictions.
 - "Demo" is displayed on MS images.
 - MS images, graphs, and data matrix tables cannot be exported.
- To use the software without these functional restrictions, or to continue using the software after 30 days, license authentication is required. When installing this software, the license management tool will be installed at the same time.
 - ▶▶ Reference "[1.2 License Authentication](#)" P.10
- When using this software for the first time after installation, you must set the location to save the data matrix temporary file. Otherwise, a data matrix cannot be created.


1.1.2 Uninstalling

NOTE Before uninstalling this software, make sure to exit the software.

1 In the [Start] menu, click [Settings] - [Apps].


2 Click [IMAGEREVEAL], then click [Uninstall].
The uninstallation window is displayed.

3 Click [Uninstall].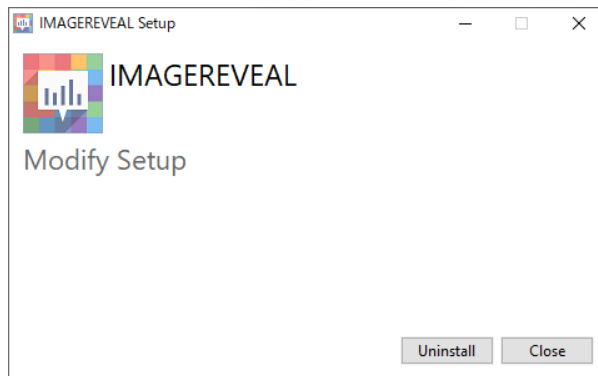

Uninstallation Window


The uninstallation starts.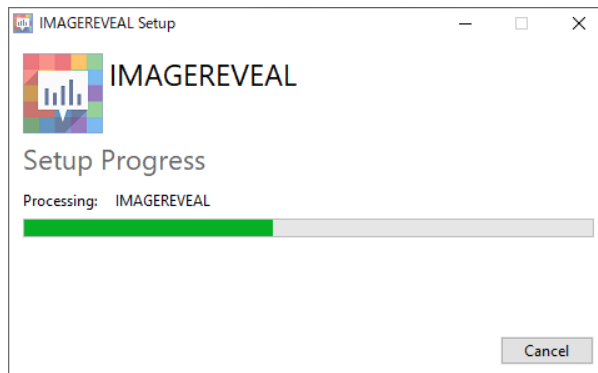

Uninstallation Progress Window

4

After the uninstallation ends successfully, click [Close] to close the uninstallation window.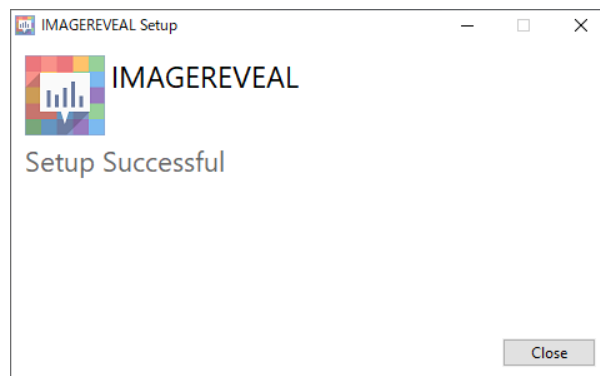

Uninstallation Complete Window

1


1.2 License Authentication

To use the MS image analysis software, license authentication is required.

You can activate or deactivate a software license by accessing the Shimadzu Software License Management System from your web browser.

NOTE The computer where this software is installed does not need to be connected to the network.

The following figure illustrates the procedure to activate a license.

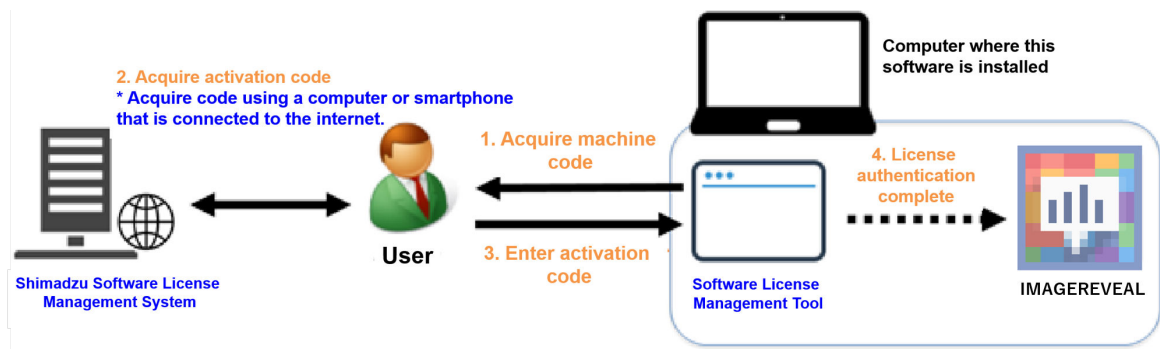License Activation Procedure

You can use the same procedure to deactivate a license.

Using the Shimadzu Software License Management System enables you to see the current usage of licenses you own. It helps you keep track of your licenses as well as manage licenses by organization and optimize the number of licenses according to the changing numbers of devices and users.

To activate or deactivate your licenses, you must register an account on the Shimadzu Software License Management System.

Register your account first by following the procedure below before using the Shimadzu Software License Management System.

1.2.1 Registering a User Account

To activate or deactivate a license, you must log in to the Shimadzu Software License Management System. Create an account for the Shimadzu Software License Management System using the following procedure.

A license certificate is also required to create an account. Have your license certificate in hand.

- NOTE
- You must register an e-mail address to create an account. Appoint a license administrator in your department and create an account for the administrator to collectively manage licenses in your department.
 - Authenticated licenses can be transferred to different accounts within the same company. You can also create a temporary account from which you can consolidate licenses to the administrator after authenticating them.
 - If you have difficulty appointing a license administrator, create an account using a group e-mail address. Even in this case, register the name of a representative.

1

Launch your web browser and access the Shimadzu Software License Management System (<https://activate.an.shimadzu.co.jp>).

2

On the login window, click [create account].

E-mail address

Password

[Forgot your password?](#)

[Log in](#)

If you do not have an account, [create account](#) from here.

Login Window

3

Enter [E-mail address], [Confirm e-mail address], and the character string from [Authentication image].

01 ▶ 02 ▶ 03 ▶ 04 ▶ 05
Input e-mail address Send e-mail Input user info Confirm contents Complete

After confirming and agreeing to the licensing and privacy policy, and consent to exchange and transfer personal data, input your e-mail address to proceed to creating an account.

E-mail address • sample@shimadzu.co.jp
Confirm e-mail address • sample@shimadzu.co.jp
Authentication image • Reload image

E-mail Input Window

4

Click [licensing and privacy policy, and consent to exchange and transfer personal data] and confirm the details.
Check [Agree to the licensing and privacy policy, and consent to exchange and transfer personal data] and click [Send].

✕ The check box can be turned on if you confirm the [licensing and privacy policy, and consent to exchange and transfer personal data](#).

Agree to the licensing and privacy policy, and consent to exchange and transfer personal data.

Send

Licensing and Privacy Policy

5

An e-mail with the subject, "[Shimadzu Corporation] Temporary registration received", is sent to the e-mail address entered in Step 2. Open your mail application and access the URL in the e-mail.

6

Enter the name, password, password confirmation, country/region, company name, and department name of the user using the system, and serial number located on the license certificate, and click [Confirm].

Input the user information.

Name *	First name :	<input type="text" value="Taro"/>	Last name :	<input type="text" value="Shimadzu"/>
E-mail address	<input type="text" value="sample@shimadzu.co.jp"/>			
Password *	<input type="password"/>			
	<small>8 or more single-byte characters</small>			
Password confirmation *	<input type="password"/>			
Country/Region *	<input type="text" value=""/>			
Company *	<input type="text"/>			
	<small>Input the company name on your business card. E.g., Shimadzu Corporation</small>			
Department	<input type="text"/>			
	<small>Input the department name on your business card. E.g., A Division, Engineering Department</small>			
	<small>To complete the account registration you must input the serial number. Input the serial number on the license certificate.</small>			
Serial number *	<input type="text"/>			

Registration Window

7

If there is no problem with the registration details on the confirmation window, click [Register].

1.2.2 Activating

If registering a license on a new computer or adding a license on a computer that has already completed license authentication, perform the following activation procedure. A license certificate is also required to activate a license. Have your license certificate in hand.

1

Power on the computer on which the IMAGEREVEAL is installed.

2


Select [Software License Management Tool], which you can find by selecting the "L" list in the Start menu and then the [License Management Tools] folder.

NOTE If the [User Account Control] window is displayed, click [Yes].

3

Click the [Acquire] button on the Activate window to acquire the machine code.

 **Hint** You can also save the machine code to a text file by clicking [Save to file].

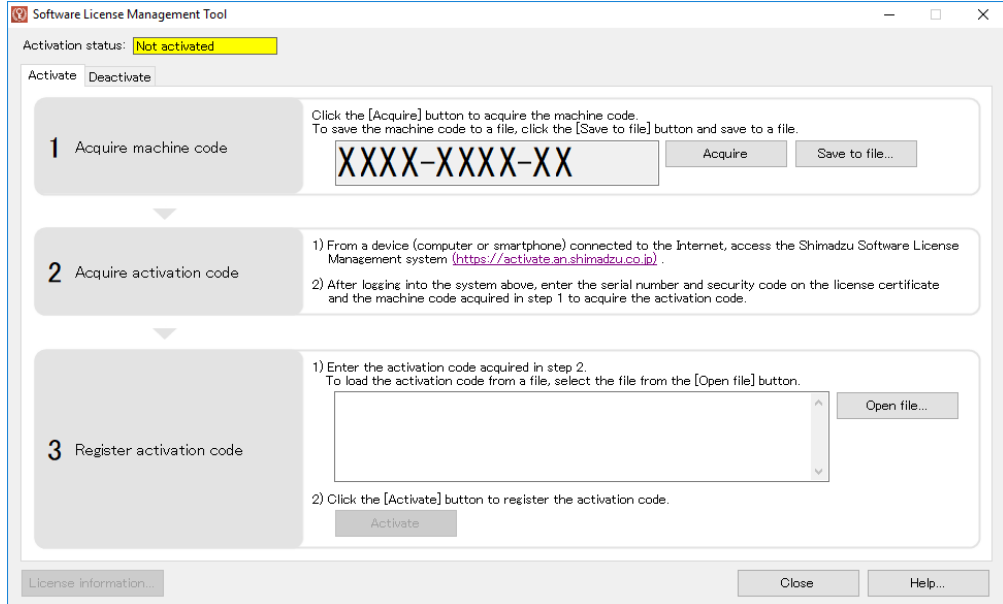Acquire Machine Code

4

From a device (computer, smartphone, etc.) connected to the Internet, access the Shimadzu Software License Management System (<https://activate.an.shimadzu.co.jp>), enter your e-mail address and password, and click [Log in].

5

Click [Activate].

Activate

Activate

6

Enter the machine code acquired in step 3 and the serial number and security code on the license certificate included with the product.

- NOTE**
- You can also enter a memo to identify the computer that is authenticating the license in the PC comment field. After entering, click [Confirm].
 - You can enter the serial number and security code of multiple licenses.

Activate

First input [machine code](#) acquired by Shimadzu Software License Management tool, and [serial number and security code](#) on the license certificate. Next click [Confirm].
* More than one license can be inputted.

Machine code * XXXX-XXXX-XX Open file

PC comment ImageExcavator

Serial number *	Security code *
L111111111	XXX

Confirm

English

© 2018 SHIMADZU Corporation. All rights reserved. [Licensing and privacy policy](#)

Activation Input Window

7

Click [Acquire] to display the activation code.

- NOTE** The activation code can also be saved to a text file by clicking [Save to file].

Activate

The activation code has been acquired.
Register the activation code on the Software License Management tool.


NNNN-NNNN-NNNN-N

Save to file

Register Activation Code

8

Enter the activation code acquired in step 7 on the Software License Management tool and click [Activate].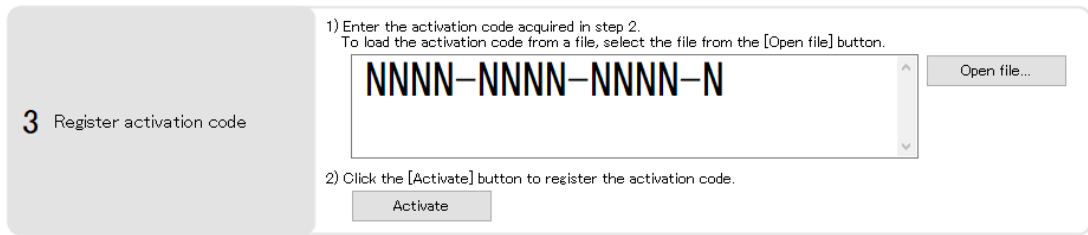

Register Activation Code

When the "Software license activation successful." message is displayed, license authentication is complete.

NOTE There is no need to restart your computer after completing license authentication.

1.2.3 Deactivating


1 Power on the computer on which the IMAGEREVEAL is installed.

2 Select [Software License Management Tool], which you can find by selecting the "L" list in the Start menu and then the [License Management Tools] folder.

NOTE If the [User Account Control] window is displayed, click [Yes].

3 Click [Deactivate] on the [Deactivate] window to acquire the deactivation code.

Hint You can also save the deactivation code to a text file by clicking [Save to file].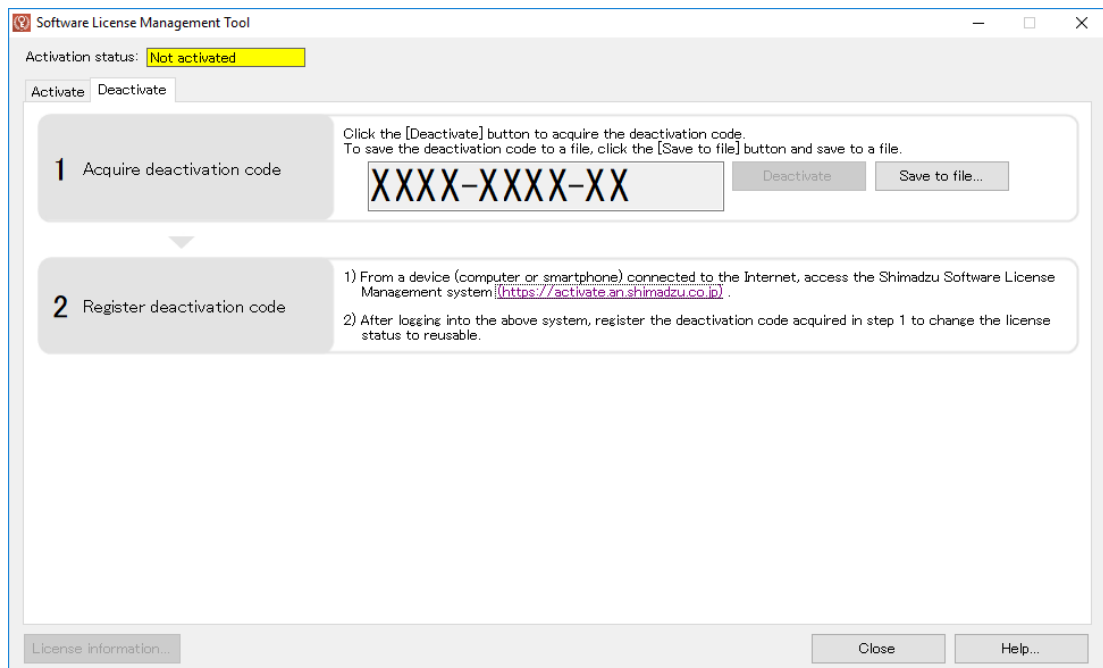

Acquire Deactivation Code

4 From a device (computer, smartphone, etc.) connected to the Internet, access the Shimadzu Software License Management System (<https://activate.an.shimadzu.co.jp>), enter your e-mail address and password, and click [Log in].

5

Click [Deactivate].


Deactivate

Deactivate

6

Enter the deactivation code acquired in step 3 and click [Register].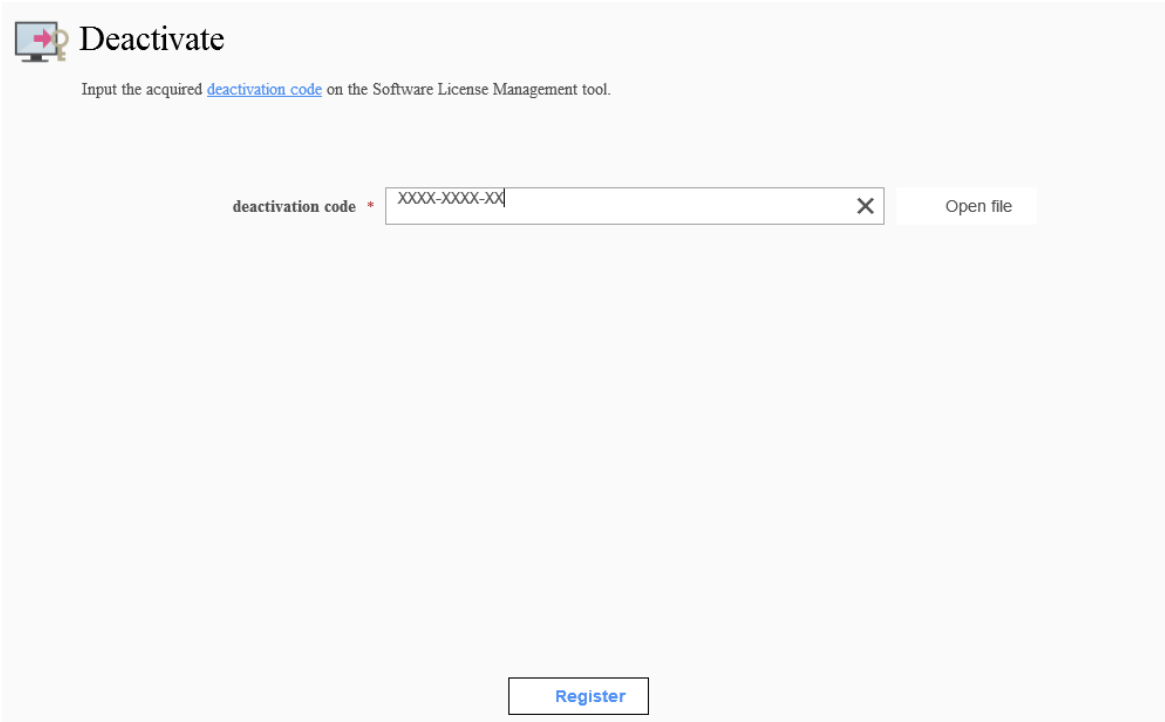

Deactivate

Input the acquired [deactivation code](#) on the Software License Management tool.

deactivation code * X Open file

Register

Register Deactivation Code

When the "The following license is reusable." message is displayed, the canceling of the license authentication is complete.

If transferring a license to another computer, activate the license on the computer receiving the transfer.

- NOTE**
- The license managed with the account logged in above step 4 is deactivated. If the license managed with another user account on the computer is activated, deactivate with another user account from above step 3. If step 3 and subsequent steps cannot be performed using another user's account, contact Shimadzu's certified service provider or your Shimadzu representative.
 - To transfer an authenticated license to another computer, you must first deactivate the license. If reusing a license on another computer, activate the license on the computer receiving the transfer after deactivating the license.
 - One license can be used on only one computer. If using multiple computers, prepare licenses for the amount of computers.
 - For troubleshooting information and frequently asked questions regarding license authentication, refer to [Frequently asked questions (FAQ)] at the top of the Shimadzu Software License Management System (<https://activate.an.shimadzu.co.jp>) window. [FAQ] is based on the installation of LabSolutions, but the contents are the same for the installation of IMAGEREVEAL. Replace "LabSolutions" with "IMAGEREVEAL" as you refer to the FAQ. If the problem persists, contact Shimadzu's certified service provider or your Shimadzu representative.

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