

# Analytical method development for widely targeted perfluoroalkyl acids (PFAAs) and their precursors in plasma using multi-gradient eluent system by LC-MS/MS

**ASMS 2017** WP-131

Jun Watanabe<sup>1</sup>, Toshikazu Minohata<sup>1</sup>, Tairo Ogura<sup>2</sup>

<sup>1</sup>Shimadzu Corporation, Japan,

<sup>2</sup>Shimadzu Scientific Instruments, Inc. Columbia, U.S.A.

# Analytical method development for widely targeted perfluoroalkyl acids (PFAAs) and their precursors in plasma using multi-gradient eluent system by LC-MS/MS

## Introduction

As perfluoroalkyl acids (PFAAs) had been widely used because of their excellent surfactant. While the product and usage of perfluorooctane sulfonic acid (PFOS), its salts and perfluorooctane sulfonyl fluoride (PFOS-F) were restricted by the Stockholm Convention in 2009, many kinds of precursors which cause PFAAs by decomposition keep being used because they are not restricted. LC-MS/MS analysis methods for PFAAs have been

developed, where a high selectivity analysis is available, but it was very difficult to implement the simultaneous analysis of PFAAs and their precursors since the wideness of their chemical property in terms of hydrophilic / hydrophobic. This time, we developed the analytical method for widely targeted PFAAs and their precursors in plasma using multi-gradient eluent system by LC-MS/MS.

## Materials and Methods

A standard mixture of PFCAs (C4-18), PFASs (C4-10), FOSAs, FTSs and PAPs, total 28 compounds, were diluted to working concentrations in methanol. Plasma sample was pretreated by WAX cartridge (final solvent: 0.05 % NH<sub>3</sub>, 0.15 % HCOOH, 80 % MeOH). Separation was achieved using Oasis WAX column (2 mmID. x 20 mL) and Triart C18 column (2 mmID. x 100 mL) maintained at 40 °C on

a UHPLC system (Shimadzu corporation, Kyoto, Japan), configured by four pumps and two 6-port valves. 500 uL of sample was inserted 250 uL water and injected. Data acquisition was performed on triple quadrupole mass spectrometer LCMS-8060 (Shimadzu Corporation, Kyoto, Japan). All samples were analyzed by multiple reaction monitoring (MRM).

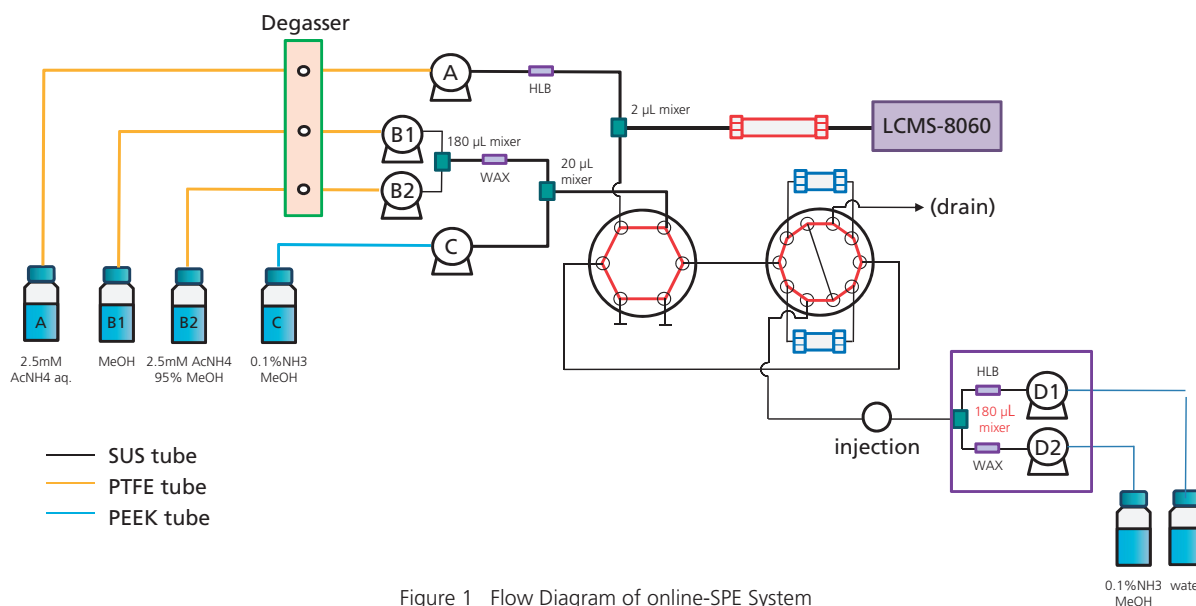


Figure 1 Flow Diagram of online-SPE System

# Analytical method development for widely targeted perfluoroalkyl acids (PFAAs) and their precursors in plasma using multi-gradient eluent system by LC-MS/MS

---

## HPLC conditions

---

Analytical Column	: YMC Triart C18 (100 mm L x 2.1 mm I.D., 3 µm)
Trap Column	: Oasis WAX (20 mm L x 2 mm I.D.)
Mobile Phase	: A: 2.5mM Ammonium Acetate in Water (Scrubber Column: OASIS HLB, 20 mm L x 2.1 mm I.D.) : B: 2.5mM Ammonium Acetate in Methanol (Scrubber Column: OASIS WAX, 20 mm L x 2.1 mm I.D.) : C: 0.1%NH <sub>4</sub> OH in MeOH
Column temperature	: 40 °C
Injection vol.	: 500 µL

---

## LCMS conditions


---

Ionization	: heated ESI negative
Nebulizing Gas Flow	: 3 L / min
Drying Gas Pressure	: 10 L / min
Heating gas flow	: 10 L / min
DL Temperature	: 150 °C
BH Temperature	: 400 °C
Interface Temperature	: 250 °C

---



# Analytical method development for widely targeted perfluoroalkyl acids (PFAAs) and their precursors in plasma using multi-gradient eluent system by LC-MS/MS

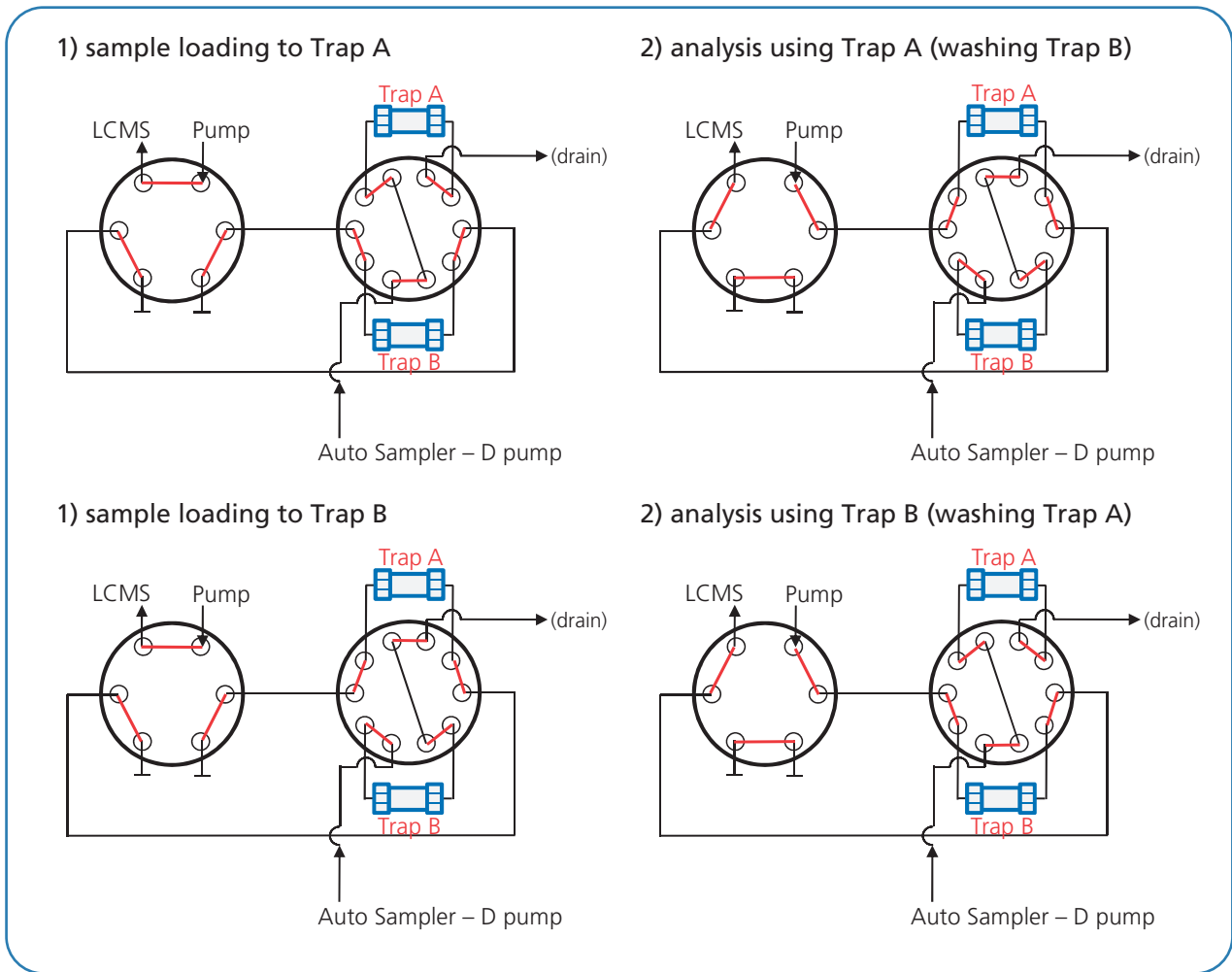



# Analytical method development for widely targeted perfluoroalkyl acids (PFAAs) and their precursors in plasma using multi-gradient eluent system by LC-MS/MS

Table 1 Calibration Range and LOD of 28 PFAAs

Compounds Name	Abbrev.	Range (pg) On column	System Blank CAL0 (pg)	LOD (ng/mL)
Perfluorobutanoic acid	PFBA	1.5-600	1	0.12
Perfluoropentanoic acid	PFPA	1.5-600	0.75	0.09
Perfluorohexanoic acid	PFHxA	1.5-600	0.75	0.09
Perfluoroheptanoic acid	PFHpA	1.5-600	0.6	0.072
Perfluorooctanoic acid	PFOA	1.5-600	0.6	0.072
Perfluorononanoic acid	PFNA	1.5-600	0.38	0.045
Perfluorodecanoic acid	PFDA	1.5-600	0.38	0.045
Perfluoroundecanoic acid	PFUnA	1.5-600	0.5	0.06
Perfluorododecanoic acid	PFDoA	1.5-600	0.38	0.045
Perfluorotridecanoic acid	PFTTrDA	1.5-600	0.38	0.045
Perfluorotetradecanoic acid	PFTeDA	1.5-600	0.5	0.06
Perfluorohexadecanoic acid	PFHxDA	1.5-600	0.75	0.09
Perfluorooctadecanoic acid	PFODA	1.5-600	0.6	0.072
Perfluorobutane sulfonate	PFBS	1.5-600	None	<0.04
Perfluorohexane sulfonate	PFHxS	1.5-600	0.3	0.036
Perfluoroheptane sulfonate	PFHpS	1.5-600	None	<0.04
Perfluorooctane sulfonate	PFOS	1.5-600	None	<0.04
Perfluorodecane sulfonate	PFDS	1.5-600	None	<0.04
N-Methyl perfluorooctane sulfonamidoacetic acid	MeFOSAA	1.5-600	None	<0.04
N-Ethyl perfluorooctane sulfonamidoacetic acid	EtFOSAA	1.5-600	None	<0.04
N-Methyl perfluorooctane sulfonamide	MeFOSA	1.5-600	None	<0.04
N-Ethyl perfluorooctane sulfonamide	EtFOSA	1.5-600	None	<0.04
4:2 Fluorotelomer sulfonic acid	4:2FTS	1.5-150	None	<0.04
6:2 Fluorotelomer sulfonic acid	6:2FTS	1.5-150	None	<0.04
8:2 Fluorotelomer sulfonic acid	8:2FTS	1.5-150	None	<0.04
6:2 Polyfluoroalkyl phosphoric acid diester	6:2diPAP	1.5-600	None	<0.04
8:2 Polyfluoroalkyl phosphoric acid diester	8:2diPAP	1.5-600	None	<0.04
Sodium bis-[2-(N-ethylperfluorooctane-1-sulfonamide)ethyl] phosphate	diSAM-PAP	1.5-600	None	<0.04

# Analytical method development for widely targeted perfluoroalkyl acids (PFAAs) and their precursors in plasma using multi-gradient eluent system by LC-MS/MS

[method A]

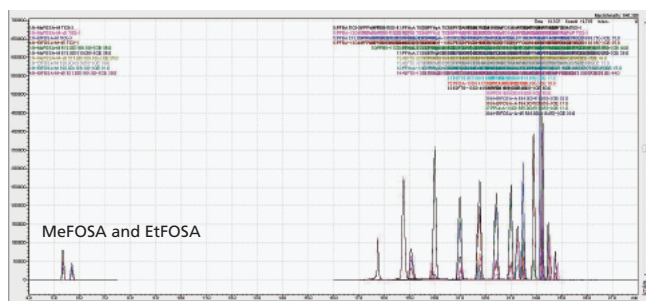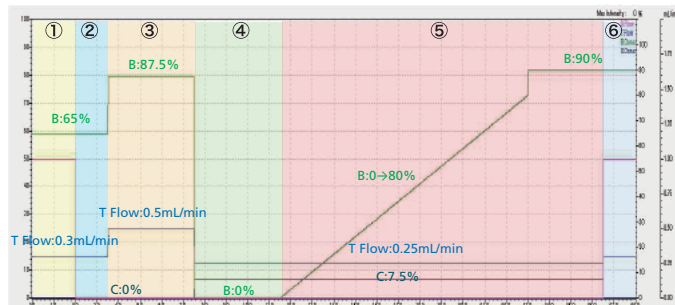

[method B]

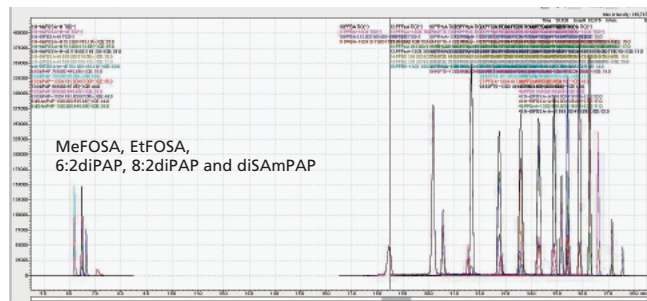
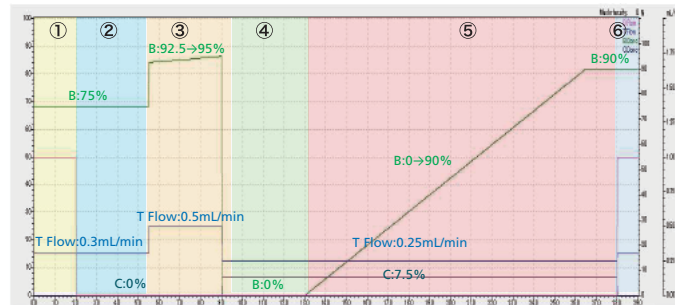


Figure 2 Gradient Program of "method A" and "method B", and MRM Chromatograms of 28 PFAAs

## Results

All of compounds were successfully ionized by negative electrospray ionization (ESI) mode. Samples were injected and loaded on trap column (WAX) with 65 % 2.5 mM NH<sub>4</sub>Ac 95 % MeOH. MeFOSA and EtFOSA were hardly retained by this condition, so these two compounds were eluted through trap column and analysis column (ODS) to mass spectrometer. After these two compounds were detected, other 26 compounds were

eluted by tertiary gradient; A: 2.5 mM NH<sub>4</sub>Ac aq., B: 2.5 mM NH<sub>4</sub>Ac MeOH, C: 0.1 % NH<sub>3</sub> MeOH, B 0 % C 7.5 % to B 80 % C 7.5%. Analysis time was set less than 30 minutes and all compounds were eluted with excellent separation. Plasma matrix samples were tested and some compounds detected with ion suppression or ion enhancement, but this method resulted good repeatability.

# Analytical method development for widely targeted perfluoroalkyl acids (PFAAs) and their precursors in plasma using multi-gradient eluent system by LC-MS/MS

Table 2 IS Recovery of 28 PFAAs comparing "method A" and "method B"

Compounds	IS recovery (%)		Compounds	IS recovery (%)	
	Method A	Method B		Method A	Method B
PFBA	58.2	35.6	PFBS	96.0	93.9
PFPA	83.1	77.1	PFHxS	73.7	62.0
PFHxA	85.2	79.6	PFHpS	92.3	86.7
PFHpA	84.5	76.6	PFOS	92.3	86.7
PFOA	96.8	88.6	PFDS	92.3	86.7
PFNA	86.4	78.6	MeFOSAA	71.6	57.2
PFDA	94.1	77.7	EtFOSAA	81.1	65.3
PFUnA	89.5	78.0	MeFOSA	82.7	76.3
PFDoA	82.3	71.4	EtFOSA	70.9	69.4
PFTTrDA	44.6	66.5	4:2FTS	44.7	23.9
PFTeDA	44.6	66.5	6:2FTS	173.6	132.4
PFHxDA	46.8	50.0	8:2FTS	142.6	80.3
PFODA	46.8	50.0	6:2diPAP	78.7	72.0
			8:2diPAP	71.8	159.9
			diSAmPAP	71.8	159.9

## Conclusions

We developed the analytical method for widely targeted PFAAs and their precursors in plasma using multi-gradient eluent system by LC-MS/MS.

## Acknowledgement

We thank Dr. Shoji F. Nakayama and Mitsuha Yoshikane (National Institute for Environmental Studies, Japan) for his help and valuable suggestions.

The product and applications in this presentation are intended for Research Use Only (RUO). Not for use in diagnostic procedures. Not available in China.

First Edition: June, 2017