

Transfer of USP normal phase method for the analysis of cortisone acetate to the Nexera UC SFC system

HPLC 2015 PSA-SFC-04

Tanaka, Kenichiro¹; Nishimura, Masayuki¹; Bollig, Brigitte²; Schad, Gesa J.²

¹ Shimadzu Scientific Instruments, Columbia, USA

PO-CON1572E

² Shimadzu Europa GmbH, Duisburg, Germany

Transfer of USP normal phase method for the analysis of cortisone acetate to the Nexera UC SFC system

Introduction

The United States Pharmacopeia (USP) monographs are widely referenced to ensure the quality of drug substances. The USP monographs offer reliable and robust methods. However, some of these methods, especially those using normal phase HPLC, require the use of harmful, even toxic solvents. Naturally, laboratories are looking to reduce the use of such solvents for health and safety, as well as cost reasons. This application note shows an example of the analysis of cortisone acetate tablets according to the USP monograph. The original normal phase method was transferred to SFC with a conventional HPLC column, Shim-pack CLC-SIL, and also a UHPLC phase, Shim-pack XR-SIL using the Nexera UC SFC system.



Figure 1: Nexera UC SFC-UV system

Method transfer to SFC

Table 1: Analytical conditions of the USP normal phase HPLC method

System Column	: Prominence : Shim-pack CLC-SIL (250 mm x 4.6 mm, 5 µm)
Mobile phase	: N-butyl chloride / water-saturated n-butyl chloride/
	Tetrahydrofuran/methanol/glacial acetic acid 95:95:14:7:6
Flow rate	: 1.0 ml/min
Column temperature	: Room temperature
Injection volume	: 15 µl
Detection	: UV 254 nm
Flow cell	: High-pressure cell for SFC

Transfer of USP normal phase method for the analysis of cortisone acetate to the Nexera UC SFC system

The USP monograph for cortisone acetate tablet analysis specifies a 250 mm x 4.6 mm silica column (L3) and n-butyl chloride/water-saturated n-butyl chloride/tetrahydrofuran /methanol/glacial acetic acid = 95/95/14/7/6 mobile phase. This normal phase method was transferred to SFC method parameters. Tables 1 and 2 show the analytical conditions for USP method and SFC method, respectively. The same column (Shim-pack CLC-SIL) was used for both of the methods.

System	: Nexera UC		
Column	: Shim-pack CLC-SIL (250 mm x 4.6 mm, 5 μm)		
Mobile phase	: A: CO ₂ (85 %)		
	B: Methanol (15 %)		
Flow rate	: 3.0 ml/min		
Column temperature	: 40 °C		
Injection volume	: 2 µl		
Detection	: UV 254 nm		
Flow cell	: High-pressure cell for SFC		

Table 2: Analytical conditions for SFC method (I)

Figure 2 shows the comparison of chromatograms of system suitability solution (methylparaben 0.01 mg/mL, cortisone acetate 0.12 mg/mL, hydrocortisone acetate 0.1 mg/mL) from a) the USP method and b) the SFC method. The analysis time was shortened from 8.5 min to 3.6 min. Table 3 shows the system suitability result. The SFC

method satisfies all of the criteria for the system suitability requirement in the USP-NF. By transferring to SFC and a CO_2 /methanol mobile phase, complex mixtures of harmful organic solvents with costly waste disposal can be changed to clean and inexpensive solvents.

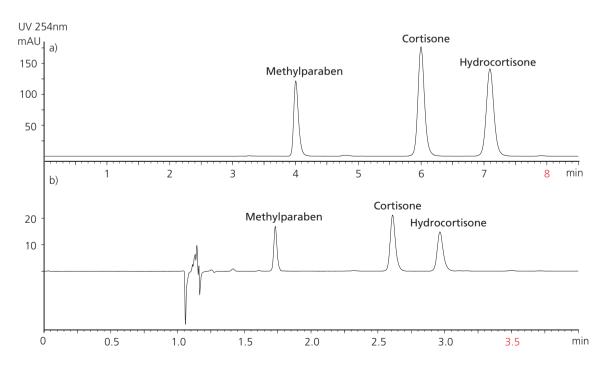


Figure 2: Chromatogram of system suitability solution with a) the USP method and b) the SFC method with conventional column

Transfer of USP normal phase method for the analysis of cortisone acetate to the Nexera UC SFC system

System suitability requirements		CLC-SIL Prominence	CLC-SIL Nexera UC
USP resolution between cortisone acetate and hydrocortisone acetate	≥ 2.2	5.33	4.64
Relative standard deviation for	≤ 2.0 %	Rt.: 0.062 %	Rt.: 0.041 %
cortisone acetate		Area: 0.133 %	Area: 0.327 %
Relative standard deviation for	≤ 2.0 %	Rt.: 0.078 %	Rt.: 0.052 %
hydrocortsione acetate		Area: 0.142 %	Area: 0.448 %

Table 3: Results of system suitability test

Method Transfer to UHPLC column

An UHPLC column with small particles was used on the Nexera UC system to further shorten the run time. Table 4 shows the analytical conditions. Figure 3 shows the comparison of chromatograms from SFC method with conventional column (Shim-pack CLC-SIL) and UHPLC column (Shim-pack XR-SIL). The runtime was decreased from 3.6 min to 0.62 min – almost 6 times shorter – while maintaining resolution. Table 5 shows the system suitability result. The SFC method with a UHPLC column satisfies all of the criteria for the system suitability requirements in the USP-NF.

Table 4: Analytical conditions for SFC method (II)

System	: Nexera UC	
Column	: Shim-pack XR-SIL (100 mm x 3.0 mm, 2.2 μm)	
Mobile phase	: A: CO ₂ (85 %)	
	B: Methanol (15 %)	
Flow rate	: 3.0 ml/min	
Column temperature	: 40 °C	
Injection volume	: 1 µl	
Detection	: UV 254 nm	
Flow cell	: High-pressure cell for SFC	

Excellence in Science

Transfer of USP normal phase method for the analysis of cortisone acetate to the Nexera UC SFC system

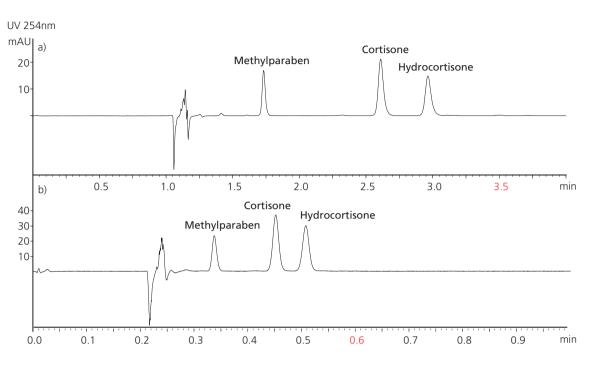


Figure3: Chromatogram of system suitability solution with a) SFC method with conventional column and b) SFC method with UHPLC column

System suitability requirements		CLC-SIL Nexera UC	CLC-SIL Nexera UC
USP resolution between cortisone acetate and hydrocortisone acetate	≥ 2.2	4.64	2.61
Relative standard deviation for	≤ 2.0 %	Rt.: 0.041 %	Rt.: 0.336 %
cortisone acetate		Area: 0.327 %	Area: 0.563 %
Relative standard deviation for	≤ 2.0 %	Rt.: 0.052 %	Rt.: 0.331 %
hydrocortsione acetate		Area: 0.448 %	Area: 0.405 %

Table 5: Results of system suitability test

Results

Figure 4 shows the comparison of analysis time and cost between the USP method and the SFC method with conventional column and UHPLC column. By transferring the original USP method to an SFC method with conventional column, analysis time and cost was shortened by a factor of 2.4 and 6.5, respectively. By transferring the original USP method to an SFC method with UHPLC column, analysis time and cost was shortened by a factor of 14 and 38, respectively. These results show that the Nexera UC enabled significant reduction of analysis time and cost without sacrificing quality of analyses.

Transfer of USP normal phase method for the analysis of cortisone acetate to the Nexera UC SFC system

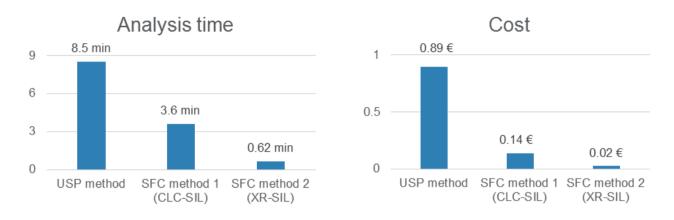


Figure 4: Comparison of time and cost using the conventional and alternative SFC method

First Edition: June, 2015



Shimadzu Corporation www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

© Shimadzu Corporation, 2015